



**KERN & Sohn GmbH**

Ziegelei 1  
D-72336 Balingen  
email: [info@kern-sohn.com](mailto:info@kern-sohn.com)

Phone: +49-[0]7433- 9933-0  
Fax: +49-[0]7433-9933-149  
Internet: [www.kern-sohn.com](http://www.kern-sohn.com)

# Installation instructions

## Tripod

### **KERN BVBP-A01**

Version 1.0  
02/2008  
GB



BVBP-A01-IA-e-0810



# KERN BVBP-A01

Version 1.0 02/2008

## Installation Instructions Tripod

Tripod with height adjustment, maximum height 1550 mm

### Scope of delivery - Assembly tripod:

	<p>1. Assemble the single parts acc. to figure</p>	<p>2. Screw-on holder of display part</p>

## Assembly tripod - Display unit:

1. Screw-off table foot of the display unit



2. Screw-off cover plate and replace it with the cover plate of the tripod



3. Screw-on cover plate



4. Screw tripod holder on the cover plate acc. to figure



5. Fix display unit on the tripod using the securing screw (1).



6. Screw-in turning knobs (2, 3) for display positioning