



KERN & Sohn GmbH

Ziegelei 1

D-72336 Balingen

E-Mail: info@kern-sohn.com

Tel: +49-[0]7433-9933-0

Fax: +49-[0]7433-9933-149

Internet: www.kern-sohn.com

Operating instruction Industrial scale

KERN FES/FEJ

Version 3.2

03/2011

GB



FES/FEJ-BA-e-1132



KERN FES/FEJ

Version 3.2 03/2011

Operating instruction Industrial scale

Contents

1	Technical data	5
2	Declaration of conformity	7
3	Basic Information (General)	8
3.1	Proper use	8
3.2	Improper Use.....	8
3.3	Warranty	8
3.4	Monitoring of Test Resources.....	9
4	Basic Safety Precautions	9
4.1	Pay attention to the instructions in the Operation Manual	9
4.2	Personnel training.....	9
5	Transport and storage	9
5.1	Testing upon acceptance	9
5.2	Packaging.....	9
6	Unpacking, Setup and Commissioning	10
6.1	Installation Site, Location of Use	10
6.2	Unpacking.....	10
6.3	Scope of delivery	11
6.4	Setup	13
6.4.1	Assembly instructions for the use of the stand (option)	15
6.5	Mains connection.....	16
6.6	Battery power supply	17
6.7	Terminal for external devices	17
6.8	Initial Commissioning.....	18
6.8.1	Power display	18
6.8.2	Bar graph display.....	19
6.8.3	Stability display.....	20
6.8.4	Balance zero display	21
6.9	Adjustment.....	21
6.9.1	Adjustment with external weight (only FES).....	21
6.9.2	Adjustment test with external weight (only FES).....	23
6.9.3	Adjustment with internal weight (only FEJ)	25
6.9.4	Adjustment test with internal weight (only FEJ)	26
6.10	GLP Log	28
6.11	Verification.....	29
6.12	Verification switch and seal	29
7	Application and configuration menu 1	31
7.1	User principle of the menu control.....	31
7.2	Menu overview	33
7.2.1	Parameter for serial interface	35

8	Configuration menu 2	37
8.1	User principle of the menu control.....	37
8.2	Menu overview	38
9	Operation	39
9.1	Keyboard overview	39
9.2	Overview of display	40
10	Weighing mode	41
10.1	Weighing.....	41
10.1.1	Taring.....	42
10.1.2	Net/gross.....	44
10.2	Weighing/parts counting.....	45
10.3	Weighing/percent determination.....	48
10.3.1	Entering the reference weight by weighing.....	48
10.3.2	Numeric entering of the reference weight.....	49
11	Adding of displayed values	52
12	Weighing with tolerance range	54
12.1	General Information.....	54
12.2	Display of the results	55
12.2.1	For 2 limits.....	55
12.2.2	For 3 or 4 limits	56
12.3	Basic settings for weighings with tolerance range.....	56
12.4	Evaluation of absolute values.....	58
12.4.1	Entering 2 limits by weighing	58
12.4.2	Entering 3 or 4 limits by weighing	61
12.4.3	Numeric entering of 2 limits	64
12.5	Evaluation with difference values	67
12.5.1	Entering 2 limits by weighing	67
12.5.2	Entering 3 or 4 limits by weighing	70
12.5.3	Numeric entering of 2 limits	70
13	Setting date and time	73
13.1	Time.....	73
13.2	Date	75
14	Description of individual functions	77
14.1	Auto sleep function	77
14.2	Auto OFF function	77
14.3	Set units function	77
14.4	Multiple range display.....	77
14.5	Interval output function	78
14.5.1	Interval setting.....	78
14.5.2	Start/Stop interval output.....	79
14.6	Input balance ID-no.	80

15	Data output	82
15.1	RS 232C interface	82
15.2	Printer interface (unidirectional data exchange).....	83
15.3	Description of interface.....	83
15.4	Data output	84
15.4.1	Format for data transmission	84
15.4.2	Signs	85
15.4.3	Data.....	85
15.4.4	Units	85
15.4.5	Result evaluation for balances with tolerance range	86
15.4.6	Data status	86
15.4.7	Interval data output	86
15.4.8	Output time.....	87
15.5	Remote control instructions	87
16	Service, maintenance, disposal	88
16.1	Cleaning	88
16.2	Service, maintenance	88
16.3	Disposal.....	88
17	Instant help	89

1 Technical data

KERN	FES 17K0.1 IPM	FES 33K0.1 IPM	FES 62K0.1 DIPM
Readability (d)	0.1g	0.1g	0.1g//1g
Verification value (e)	1g	1g	-
Weighing range (max)	17kg	33kg	6.2kg//62kg
Accuracy class	II	II	II
Warming up time (operating temperature)	2 h	2 h	2 h
Taring range (subtractive)	17kg	33kg	62kg
Minimum load (Min)	5g	5g	-
Minimum piece weight	0.1 g	0.1g	0.1g
Reference quantities	5,10, 30, 100	5,10, 30, 100	5, 10, 30, 100
Reproducibility	0,1	0,1	0.1g//1g
Linearity	± 0.3g	± 0.3 g	± 0.3g / 1g
Recommended adjusting weight, not included (class)	10kg + 5kg (F1)	20kg + 10kg (E2)	50 kg (F1)
Stabilization time	3 sec.	3 sec.	3 sec.
Weighing plate, stainless steel	350x400x170mm	350x400x170mm	350x400x170mm
Net weight (kg)	17kg	17kg	17kg
Units	g, kg, ct		
Humidity of air	max. 80 % relative (not condensing)		
Permissible ambient condition	5° C to 35° C		
Protection type	IP65 in battery operation		
Dimensions of the balance(B x D x H)	350 x 520 x 170 mm		
Vibration filter	4		
Mains connection	Mains adaptor 230 V, 50/60 Hz ; 12 V DC balance, 600 mA		
Rechargeable battery (optional)	Operating time ca. 6h. / charging time ca. 12h		

KERN	FEJ 17K0.1 IPM	FEJ 33K0.1 IPM	FEJ 62K0.1 DIPM
Readability (d)	0.1g	0.1g	0.1g//1g
Verification value (e)	1g	1g	-
Weighing range (max)	17kg	33kg	6.2kg//62kg
Accuracy class	II	II	II
Warming up time (operating temperature)	2 h	2 h	2 h
Taring range (subtractive)	17kg	33kg	62kg
Minimum load (Min)	5g	5g	-
Minimum piece weight	0.1 g	0.1g	0.1g
Reference quantities	5,10, 30, 100	5,10, 30, 100	5, 10, 30, 100
Reproducibility	0,1	0,1	0.1g//1g
Linearity	± 0.3g	± 0.3 g	± 0.3g / 1g
Adjustment weight	internal	internal	internal
Stabilization time	3 sec.	3 sec.	3 sec.
Weighing plate, stainless steel	350x400x170mm	350x400x170mm	350x400x170mm
Net weight (kg)	17kg	17kg	17kg
Units	g, kg, ct		
Humidity of air	max. 80 % relative (not condensing)		
Permissible ambient condition	5° C to 35° C		
Protection type	IP65 in battery operation		
Dimensions of the balance(B x D x H)	350 x 520 x 170 mm		
Vibration filter	4		
Mains connection	Mains adaptor 230 V, 50/60 Hz ; 12 V DC balance, 600 mA		
Rechargeable battery (optional)	Operating time ca. 6h. / charging time ca. 12h		

2 Declaration of conformity



KERN & Sohn GmbH

D-72322 Balingen-Frommern

Postfach 4052

E-Mail: info@kern-sohn.de

Tel: 0049-[0]7433- 9933-0

Fax: 0049-[0]7433-9933-149

Internet: www.kern-sohn.de

Konformitätserklärung

EC-Konformitätserklärung

EC- Déclaration de conformité

EC-Dichiarazione di conformità

EC- Declaração de conformidade

EC-Deklaracja zgodności

EC-Declaration of -Conformity

EC-Declaración de Conformidad

EC-Conformiteitverklaring

EC- Prohlášení o shode

ЕС-Заявление о соответствии

D	Konformitäts- erklärung	Wir erklären hiermit, dass das Produkt, auf das sich diese Erklärung bezieht, mit den nachstehenden Normen übereinstimmt.
GB	Declaration of conformity	We hereby declare that the product to which this declaration refers conforms with the following standards.
CZ	Prohlášení o shode	Tímto prohlašujeme, že výrobek, kterého se toto prohlášení týká, je v souladu s níže uvedenými normami.
E	Declaración de conformidad	Manifestamos en la presente que el producto al que se refiere esta declaración está de acuerdo con las normas siguientes
F	Déclaration de conformité	Nous déclarons avec cela responsabilité que le produit, auquel se rapporte la présente déclaration, est conforme aux normes citées ci-après.
I	Dichiarazione di conformità	Dichiariamo con ciò che il prodotto al quale la presente dichiarazione si riferisce è conforme alle norme di seguito citate.
NL	Conformiteit- verklaring	Wij verklaren hiermede dat het product, waarop deze verklaring betrekking heeft, met de hierna vermelde normen overeenstemt.
P	Declaração de conformidade	Declaramos por meio da presente que o produto no qual se refere esta declaração, corresponde às normas seguintes.
PL	Deklaracja zgodności	Niniejszym oświadczamy, że produkt, którego niniejsze oświadczenie dotyczy, jest zgodny z poniższymi normami.
RUS	Заявление о соответствии	Мы заявляем, что продукт, к которому относится данная декларация, соответствует перечисленным ниже нормам.

Electronic Balance: KERN PES, PEJ, FES, FEJ

Mark applied	EU Directive	Standards
	2004/108/EC	EN 55022 (2006)
	2006/95/EC	EN 60950 (2001)

Date: 07.10.2008

Signature:

**KERN & Sohn GmbH
Management**

KERN & Sohn GmbH, Ziegelei 1, D-72336 Balingen, Tel. +49-[0]7433/9933-0, Fax +49-[0]7433/9933-149

3 Basic Information (General)

3.1 Proper use

The balance you purchased is intended to determine the weighing value of material to be weighed. It is intended to be used as a “non-automatic” balance, i.e. the material to be weighed is manually and carefully placed in the centre of the weighing plate. As soon as a stable weighing value is reached the weighing value can be read.

3.2 Improper Use

Do not use balance for dynamic weighings. In the event that small quantities are removed or added to the material to be weighed, incorrect weighing results can be displayed due to the “stability compensation” in the balance. (Example: Slowly draining fluids from a container on the balance.)

Do not leave permanent load on the weighing plate. This may damage the measuring system.

Impacts and overloading exceeding the stated maximum load (max) of the balance, minus a possibly existing tare load, must be strictly avoided. This could cause damage to the balance.

Never operate balance in explosive environment. The serial version is not explosion protected. The structure of the balance may not be modified. This may lead to incorrect weighing results, safety-related faults and destruction of the balance.

The balance may only be used according to the described conditions. Other areas of use must be released by KERN in writing.



The appliance may only be opened by trained service technicians according to KERN standards.

Before opening, disconnect appliance from power supply!

Warranty claims will be voided when appliance is opened.



The **FEJ** weighing system may not be used in explosive areas or areas with explosive substances.

3.3 Warranty

Warranty claims shall be voided in case

- Our conditions in the operation manual are ignored
- The appliance is used outside the described uses
- The appliance is modified or opened
- Mechanical damage and damage caused by media, liquids
- Natural wear and tear
- The appliance is improperly set up or incorrectly electrically connected
- The measuring system is overloaded

3.4 Monitoring of Test Resources

In the framework of quality assurance the measuring-related properties of the balance and, if applicable, the testing weight, must be checked regularly. The responsible user must define a suitable interval as well as type and scope of this test. Information is available on KERN's home page (www.kern-sohn.com) with regard to the monitoring of balance test substances and the test weights required for this. In KERN's accredited DKD calibration laboratory test weights and balances may be calibrated (return to the national standard) fast and at moderate cost.

4 Basic Safety Precautions

4.1 Pay attention to the instructions in the Operation Manual

Carefully read this operation manual before setup and commissioning, even if you are already familiar with KERN balances.

4.2 Personnel training

The appliance may only be operated and maintained by trained personnel.

5 Transport and storage

5.1 Testing upon acceptance

When receiving the appliance, please check packaging immediately, and the appliance itself when unpacking for possible visible damage.

5.2 Packaging

Keep all parts of the original packaging in case you need to return the appliance. Only use original packaging for returning.

Before sending, disconnect all connected cables and loose/movable parts.

Attach possibly existing transport safeguards. Secure all parts, e.g. glass windshield, weighing plate, power unit etc., to prevent slipping and damage.

6 Unpacking, Setup and Commissioning

6.1 Installation Site, Location of Use

The balances are designed in a way that reliable weighing results are achieved in common conditions of use.

You will work accurately and fast, if you select the right location for your balance.

Therefore, observe the following for the installation site:

- Place the balance on a firm, level surface;
- Avoid extreme heat as well as temperature fluctuation caused by installing next to a radiator or in the direct sunlight;
- Protect the balance against direct draughts due to open windows and doors;
- Avoid jarring during weighing;
- Protect the balance against high humidity, vapours and dust;
- Do not expose the device to extreme dampness for longer periods of time. Non-permitted condensation (condensation of air humidity on the appliance) may occur if a cold appliance is taken to a considerably warmer environment. In this case, acclimatize the disconnected appliance for ca. 2 hours at room temperature.
- Avoid static charging of the material to be weighed, weighing container and windshield.

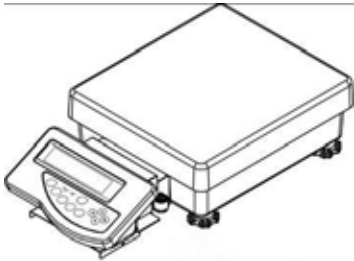
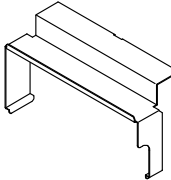
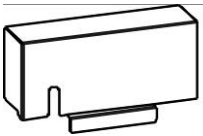
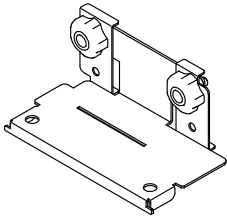
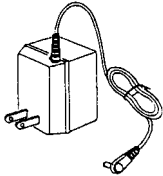
If electro-magnetic fields or static charge occur, or if the power supply is unstable major deviations on the display (incorrect weighing results) are possible. In that case, the location must be changed.

6.2 Unpacking

Carefully remove the balance from the packaging, remove plastic cover and setup balance at the intended workstation.

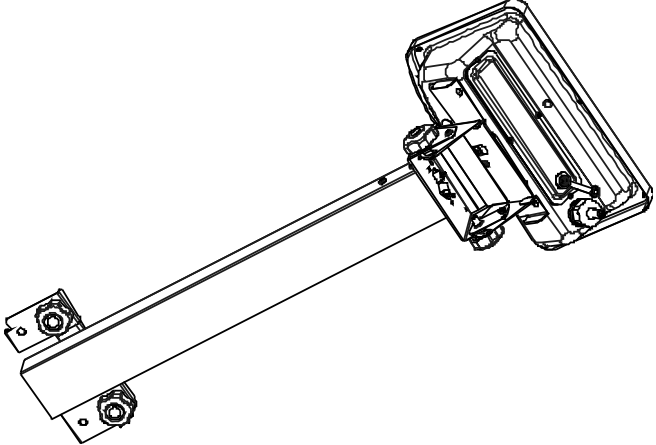
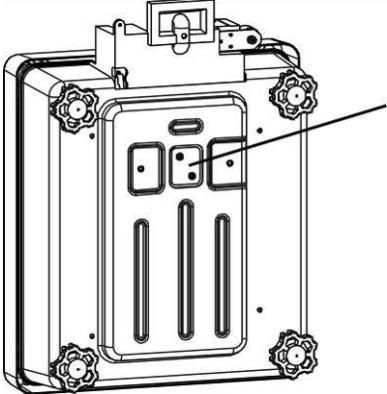
6.3 Scope of delivery

Serial accessories:

1. Balance	
2. Cable box cover	
3. Cover display holder	
4. Display holder	
5. Mains power supply	
6. Operating Manual	

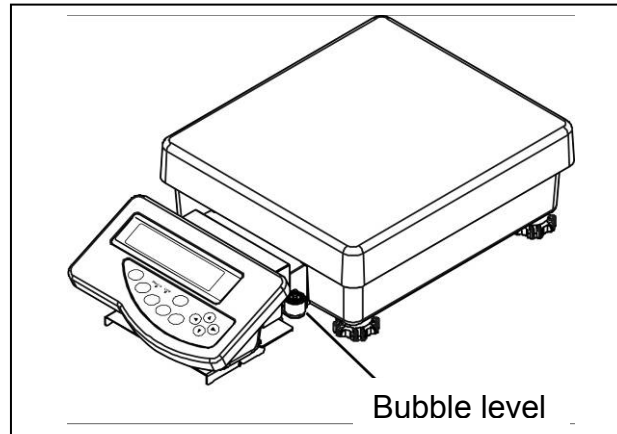
Accessories optional:

English

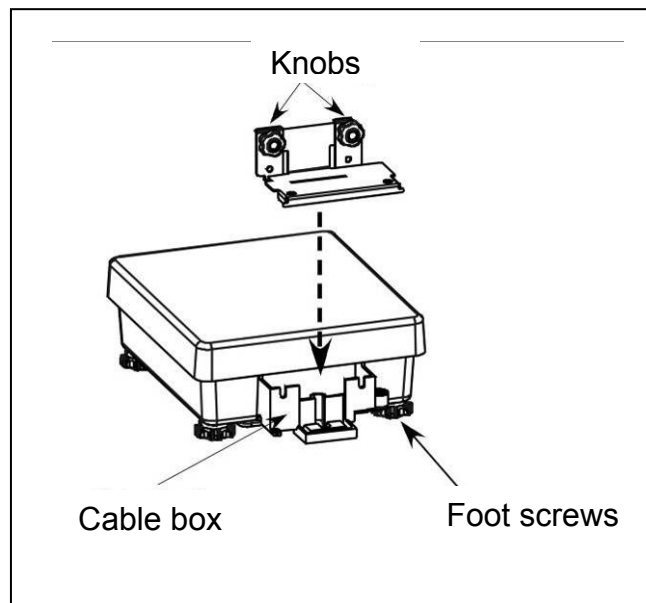
<p>1. Stand (FEJ-A05)</p>	
<p>2. Hook for underfloor weighing (FEJ-A06)</p>	 <p>Remove closing cover in order to screw-on the hook. To protect against dust and humidity put it again after the underfloor-weighing.</p>
<p>3. Relay output (FEJ-A07)</p>	

6.4 Setup

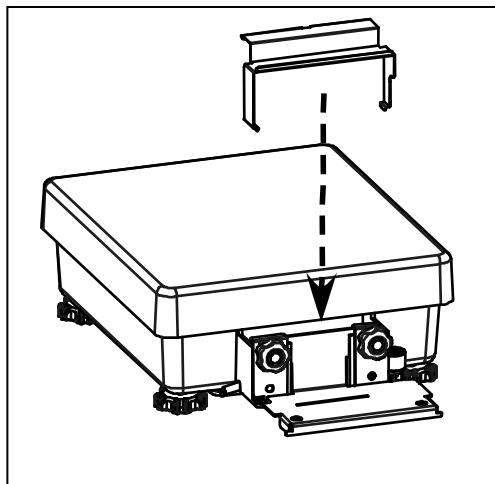
1. Level balance with foot screws until the air bubble of the water balance is in the prescribed circle.



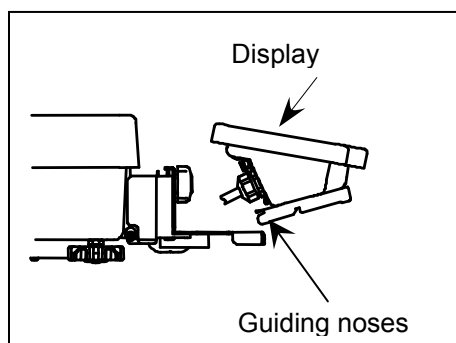
2. Remove weighing plate



3. Release the knobs on the display holder almost completely
4. Suspend the display holder on the cable box from above
5. Turn in knobs completely in order to fix them
6. Draw the cable into the cable box, bend and store it, and let 15 cm remain outside.

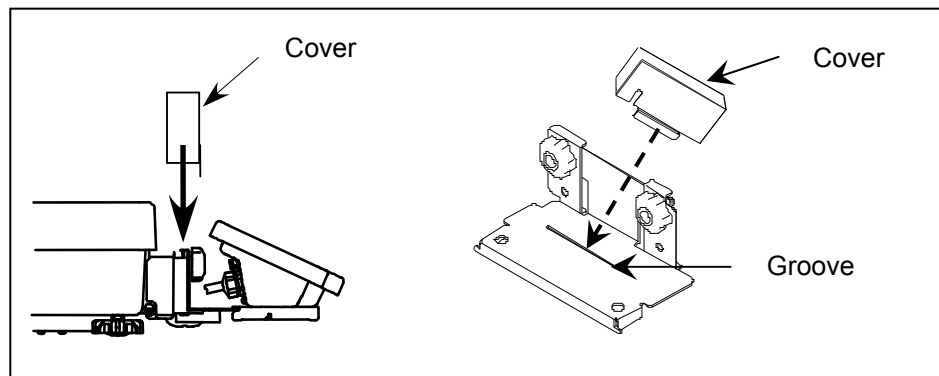


7. Apply the cable box cover from above, bending the sides slightly outwards.
8. Hold the display a little inclined upwards and slide the lateral guiding noses under the lateral grooves of the display holder.

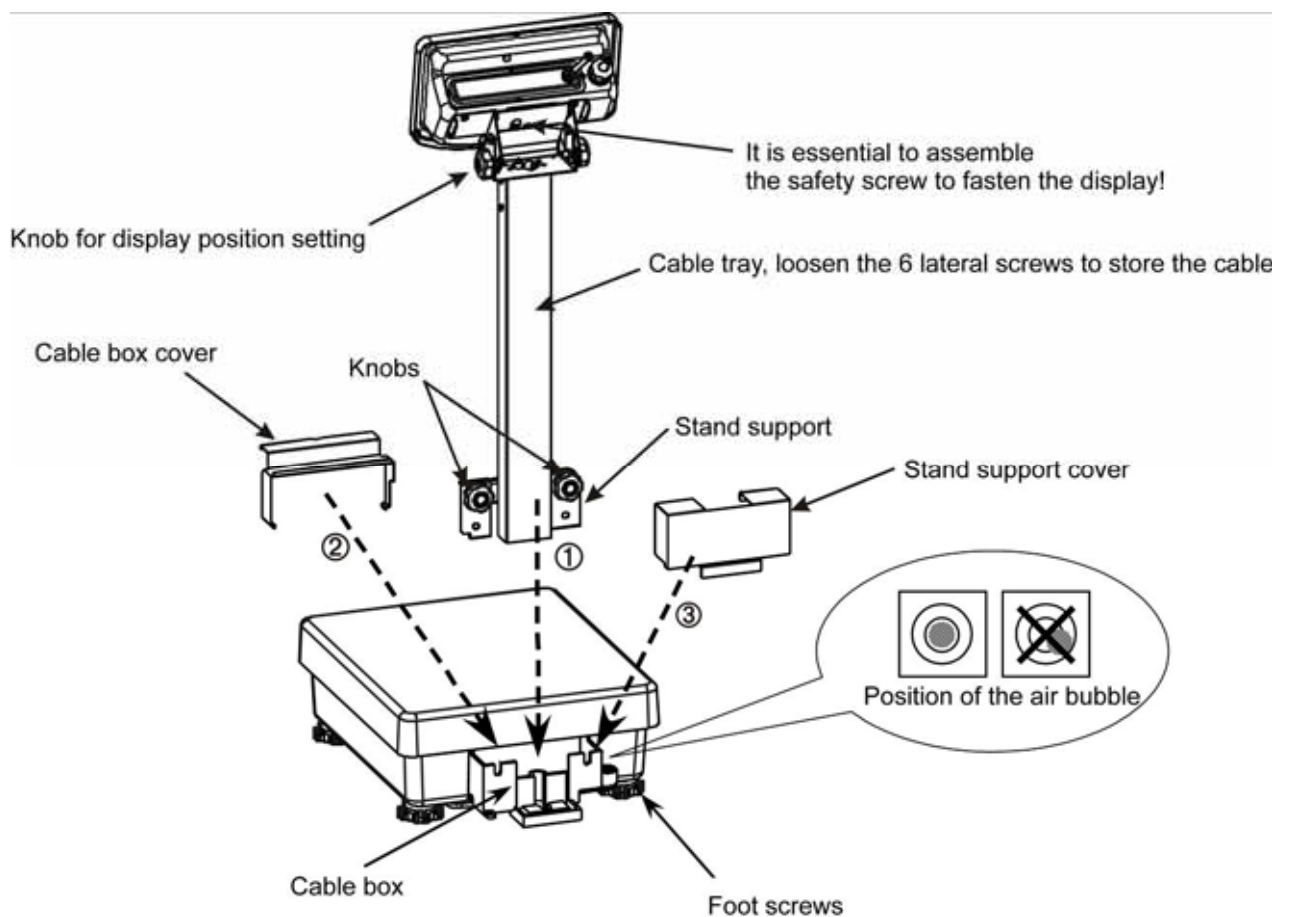


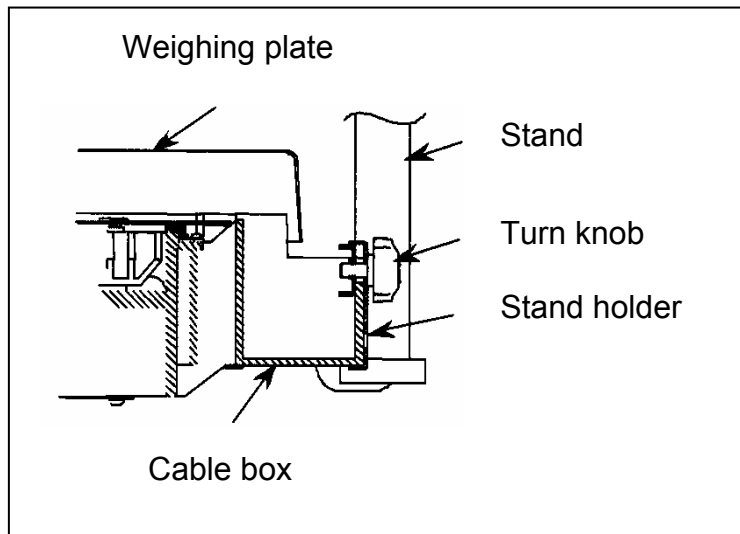
9. Re-install the weighing plate

10. Insert the display holder cover from the top above the two knobs of the display holder. It is locked in the groove on the display holder.



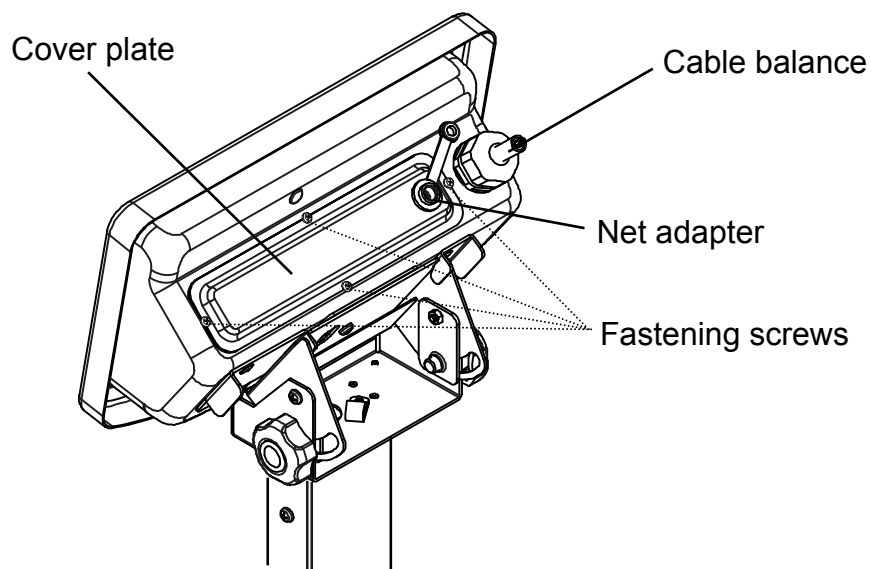
6.4.1 Assembly instructions for the use of the stand (option)



Lateral view:**6.5 Mains connection**

Power is supplied via the external mains power supply. The stated voltage value must be the same as the local voltage.

Only use original KERN mains power supplies. Using other makes requires consent by KERN.






6.6 Battery power supply

The optionally supplied battery is charged with the supplied power supply.

Before the first use, the battery should be charged by connecting it to the mains power supply for at least 12 hours. The operating time of the battery is about. 6h. Charging time until complete recharging ca. 12h.

In the menu you can activate the AUTO-OFF function [9 R.P. !]. After 3 min without load change the balance switches automatically off in order to spare the battery. Whilst in battery mode the balance is IP 65 protected.

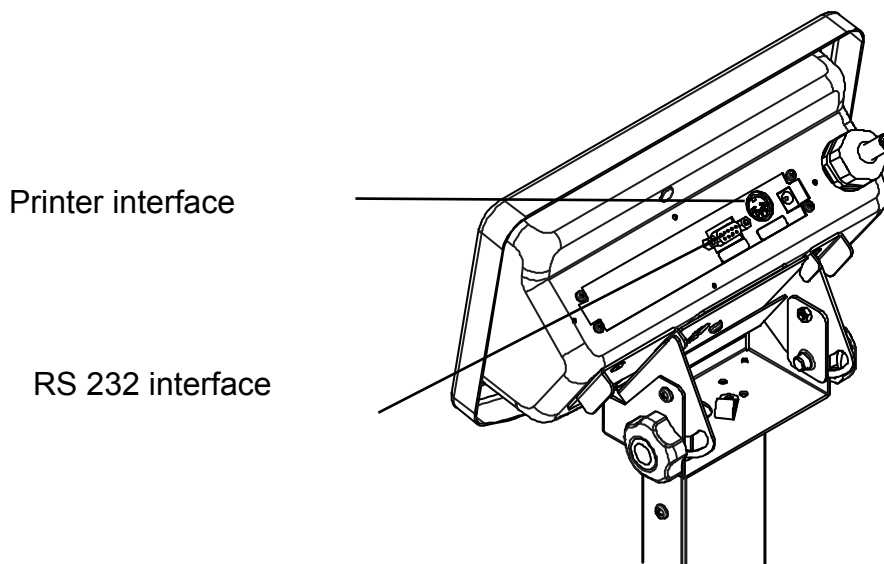
When the balance is in battery mode the following symbols appear on the display:

	Battery charge sufficient
	Battery very low. To charge the battery, connect it to the mains as soon as possible (re-calibration not possible).
	Voltage has dropped below prescribed minimum (7V). Plug in the mains adapter, to charge the balance via the electrical network (12h).

6.7 Terminal for external devices

The output for the external devices is situated under the cover plate on the rear of the display. To detach the cover plate, remove the fastening screws (see fig. chap. 6.5).

Caution: Doing this, the IP 65 protection provided in battery mode will be lost.



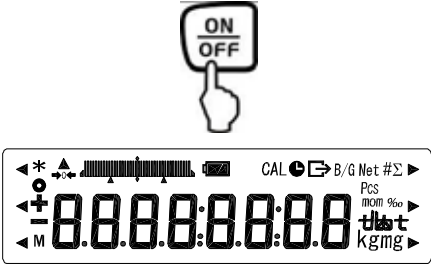


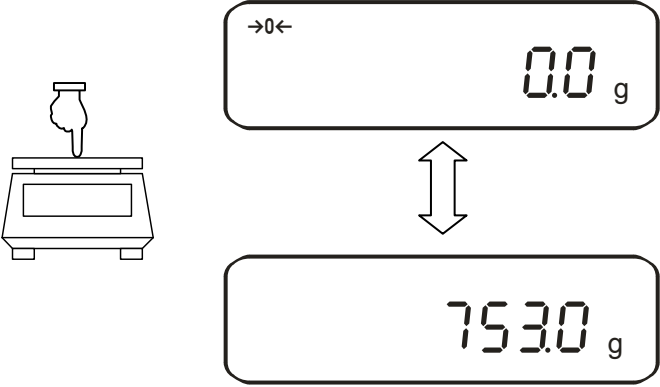

6.8 Initial Commissioning

In order to obtain exact results with the electronic balances, your balance must have reached the operating temperature (see warming up time chap. 1). During this warming up time, the balance must be connected to the power supply (mains, accumulator or battery).

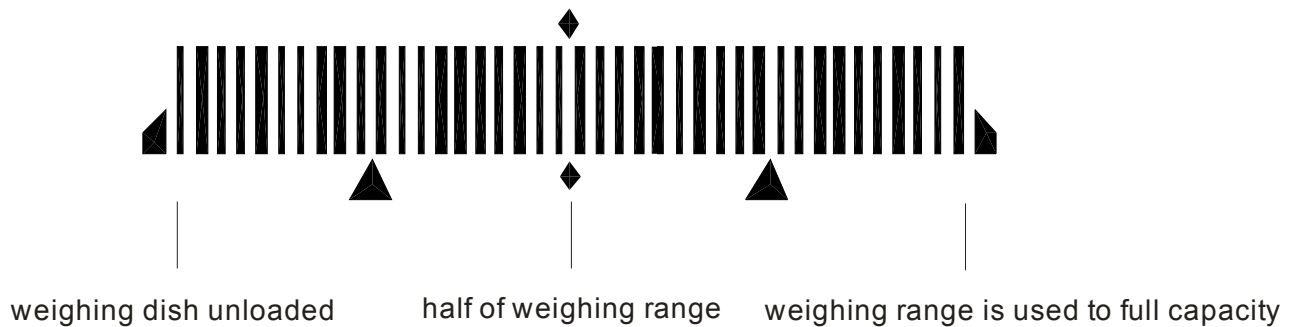
The accuracy of the balance depends on the local acceleration of gravity.

Please be sure to observe the information in the chapter on adjusting in **chap. 6.8**.

6.8.1 Power display

 <p>The balance will carry out a self-test</p>	<p>Supply balance with power via the mains power supply.</p> <p>Balance is in stand-by mode (green LED is on).</p> <p>Use the -key to switch the balance on.</p>
	<p>If there is an object on the weighing plate when the balance is switched on there is a flashing display „on 0“. In that case, remove object from the weighing plate.</p>
	<p>By pressing lightly it is possible to check whether the balance display changes.</p>
<p>Stand-by ■</p>	<p>Use the -key to switch the balance off. The balance is now in stand-by mode again (green LED is on).</p>

6.8.2 Bar graph display



The weighing range of the balance is divided into 40 graphic cuboids. Zero (0) will appear on the graphic display if there is no weighing value on the balance. 20 graphic cuboids are displayed if the balance is loaded up to one half of its weighing range.

In configuration menu 1 (chapter 7) you can activate/deactivate the bar graph display.

Parameter selection:

g b.c.	0	Hide bar graph
	* 1	Show bar graph

Bar graph display in the respective operating modes:

Bar graph display	Operating mode
	Weighing mode Tolerance weighing with 1 or 2 limits
	Tolerance weighing with 3 or 4 limits

NOTE:

If tare weighing is being carried out, the graphic weight display will continue to indicate the number of cuboids of the tare weight.

6.8.3 Stability display

Stable



Unstable





If the display shows the stability display **[o]** the balance is in a stable status. The **[o]** indication disappears if the condition is unstable.

6.8.4 Balance zero display

Environmental influences can lead to the exact figure of “000.0” not being displayed in spite of an empty weighing dish. It is, however, possible to reset your balance to zero at any time and thus ensure that weighing really does commence at zero. Setting to zero when a weight is applied is only possible within a certain type-dependent range. In the event that the balance cannot be reset to zero with an applied weight, this range has been exceeded.

[o - Err] will appear on the display.

Operation	Display
If the balance does not show exactly zero although the pan scale is unloaded, press the  -button and the balance will be reset to zero.	
After a brief waiting time your balance is reset to zero. In addition to this, the sign for the balance zero setting will be displayed [→0←].	

6.9 Adjustment





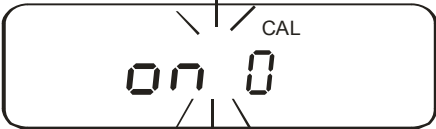




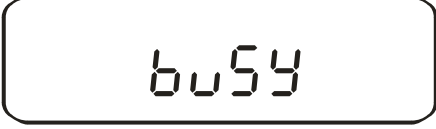
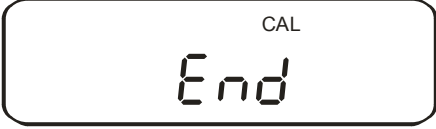

As the acceleration value due to gravity is not the same at every location on earth, each balance must be coordinated - in compliance with the underlying physical weighing principle - to the existing acceleration due to gravity at its place of location (only if the balance has not already been adjusted to the location in the factory). This adjustment process must be carried out for the first commissioning, after each change of location as well as in case of fluctuating environment temperature. To receive accurate measuring values it is also recommended to adjust the balance periodically in weighing operation.

6.9.1 Adjustment with external weight (only FES)

Procedure when adjusting:

Observe stable environment conditions. A warming-up time (see cap. 1) for stabilisation is necessary. Ensure that there are no objects on the weighing plate.

At verified balances, the adjustment is locked by a switch. In order to adjust, open the locking switch see chap.6.11.

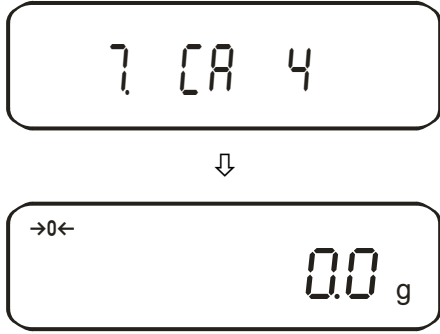
Operation	Display
Activate function [] [A. 3] (see chap. 7).	 ↓ 
 Zero point will be saved.	 ↓  If the display indicates [PuSH. F], press the  -key
Carefully place adjusting weight in the centre of the weighing plate Adjustment process is started. The process of adjustment is completed. Remove adjusting weight, balance will return into weighing mode automatically. In case of an adjustment error or incorrect adjusting weight the display will show [- Err]; repeat adjustment process.	 ↓ If the display indicates [PuSH. F], press the  -key  ↓  ↓  ↓ 


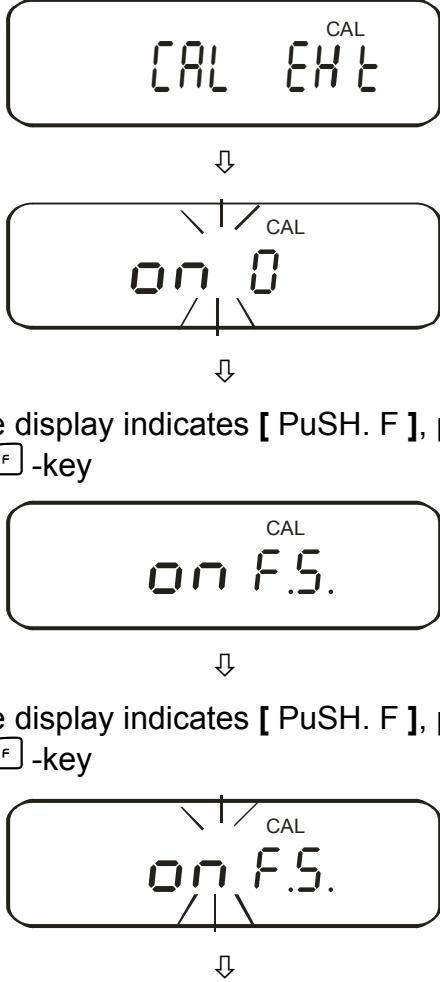
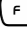

6.9.2 Adjustment test with external weight (only FES)

During adjustment tests the balance automatically compares the saved value of the adjustment weight with the actual value. This is only a check, i.e. no values are changed.

Procedure:

Observe stable environmental conditions. A warming up time (see chap. 1) is required for stabilization. Ensure that there are no objects on the weighing plate.

Operation	Display
Activate function [?] [R. 4] (see chap. 7).	

<p>Start of the adjustment test:</p>  <p>Carefully place adjusting weight in the centre of the weighing plate</p> <p>The difference between the saved value and the measured value is displayed.</p>	 <p>If the display indicates [PuSH. F], press the  -key</p> <p>If the display indicates [PuSH. F], press the  -key</p>
---	--

Take away adjustment weight.

Press any key; the adjustment process is cancelled and the balance returns to weighing mode.

dlFF



-0.02_g



CAL
End



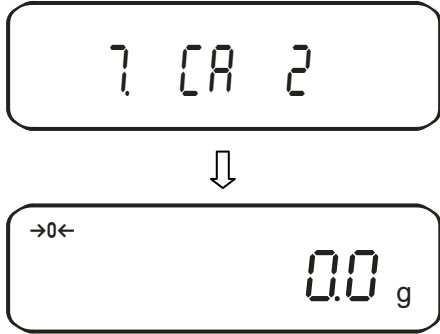

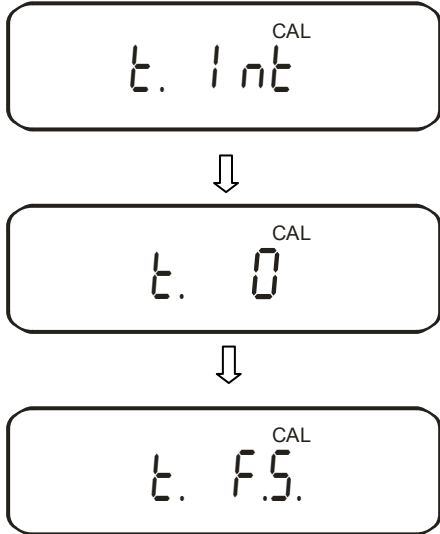
→0←
0.0_g

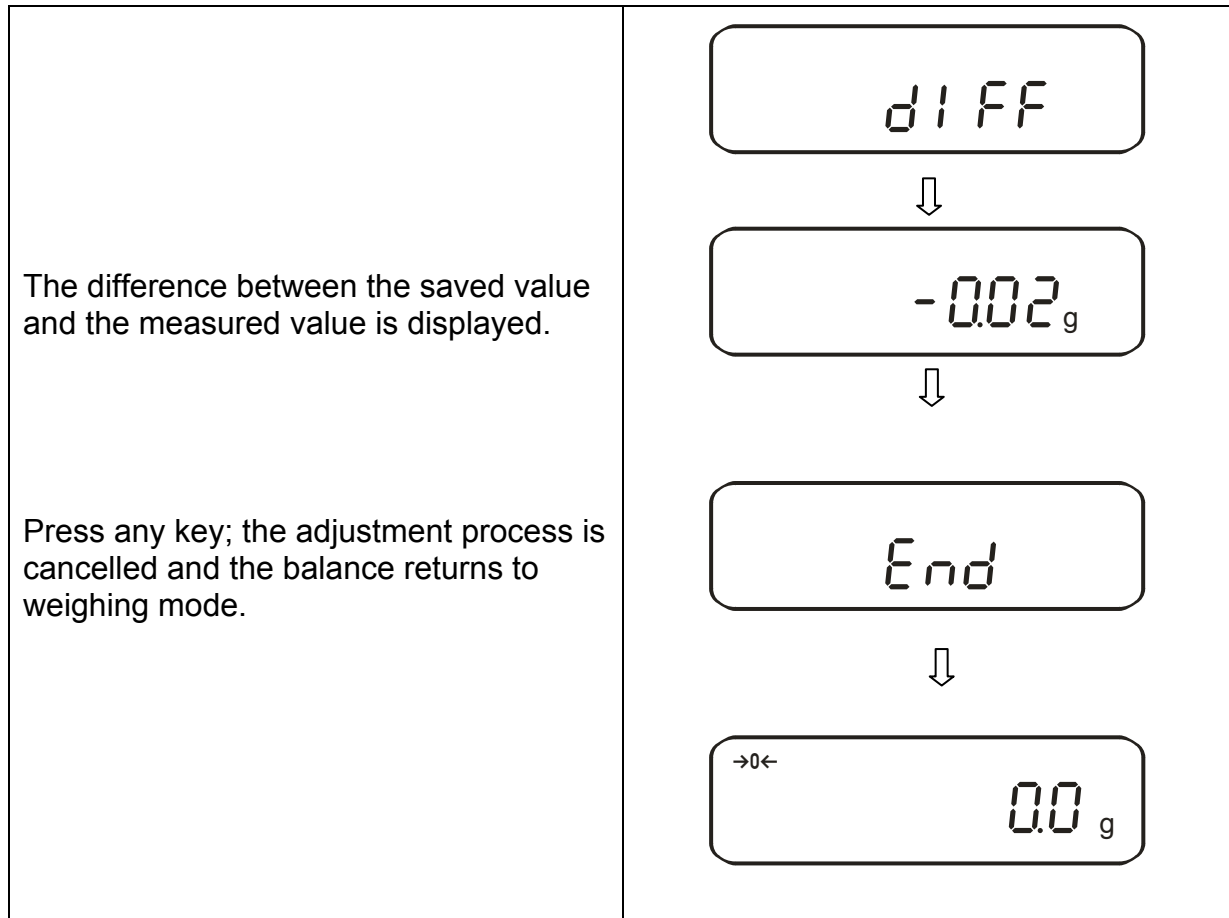
6.9.4 Adjustment test with internal weight (only FEJ)

During adjustment tests the balance automatically compares the saved value of the adjustment weight with the actual value. This is only a check, i.e. no values are changed.

Procedure:

Observe stable environmental conditions. A warming up time (see chap. 1) is required for stabilization. Ensure that there are no objects on the weighing plate.

Operation	Display
Activate function [7] [CAL. 2] (see chap. 7).	
Start of the adjustment test:  The test is carried out automatically.	



Error messages during adjustment function:

1. **3-Err:** The weighing plate is not empty -> Remove goods to be weighed from weighing plate
2. **7-Err:** Battery capacity too low -> connect balance to power adapter (load battery)

Should other error messages occur, switch balance off and then on again. If the error message remains inform manufacturer.

6.10 GLP Log

Quality assurance systems require printouts of weighing results as well as of correct adjustment of the balance stating date and time as well as balance identification. The easiest way is to have a printer connected.

Activate function (**E.GLP – 1**), see chap.7

Printout example:

DATE	04.11.2008	Date
TIME:	08:42	Time
GOTTL.KERN&SOHN		Company
TYPE	FEJ33K01IPM	Model
S/N	Dxxxxxxxx	Serial no.
ID	1234	Balance identification no.
CAL.INTERVAL		
REF:		
	033000.0 g	Adjustment weight
COMPLETE		
DATE:	04.11.2008	
TIME:	08:42	
SIGNATURE		prepared by

6.11 Verification

General introduction:

According to EU directive 90/384/EEC balances must be verified if they are used as follows (legally controlled area):

- a) For commercial transactions if the price of goods is determined by weighing
- b) For the production of medicines in pharmacies as well as for analyses in the medical and pharmaceutical laboratory
- c) For official purposes.
- d) For manufacturing final packages.

In cases of doubt, please contact your local trade in standard.

Verification instructions

An EU type approval exists for balances described in their technical data as verifiable. If a balance is used where obligation to verify exists as described above, it must be verified and re-verified in regular intervals.

Re-verification of a balance is carried out according to the respective national regulations. The validity for verification of balances in Germany is e.g. 2 years.

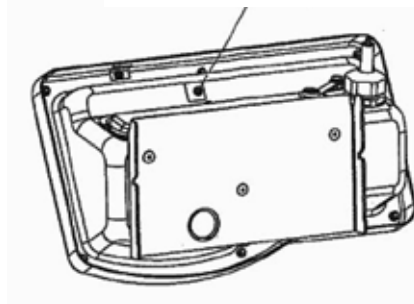
The legal regulation of the country where the balance is used must be observed!

6.12 Verification switch and seal

The calibration switch is located on the back of the display. Access to the calibration switch by removing the seal and the rubber plug (see drawing). Prior to calibrating, the calibration switch must be moved into the calibration position.

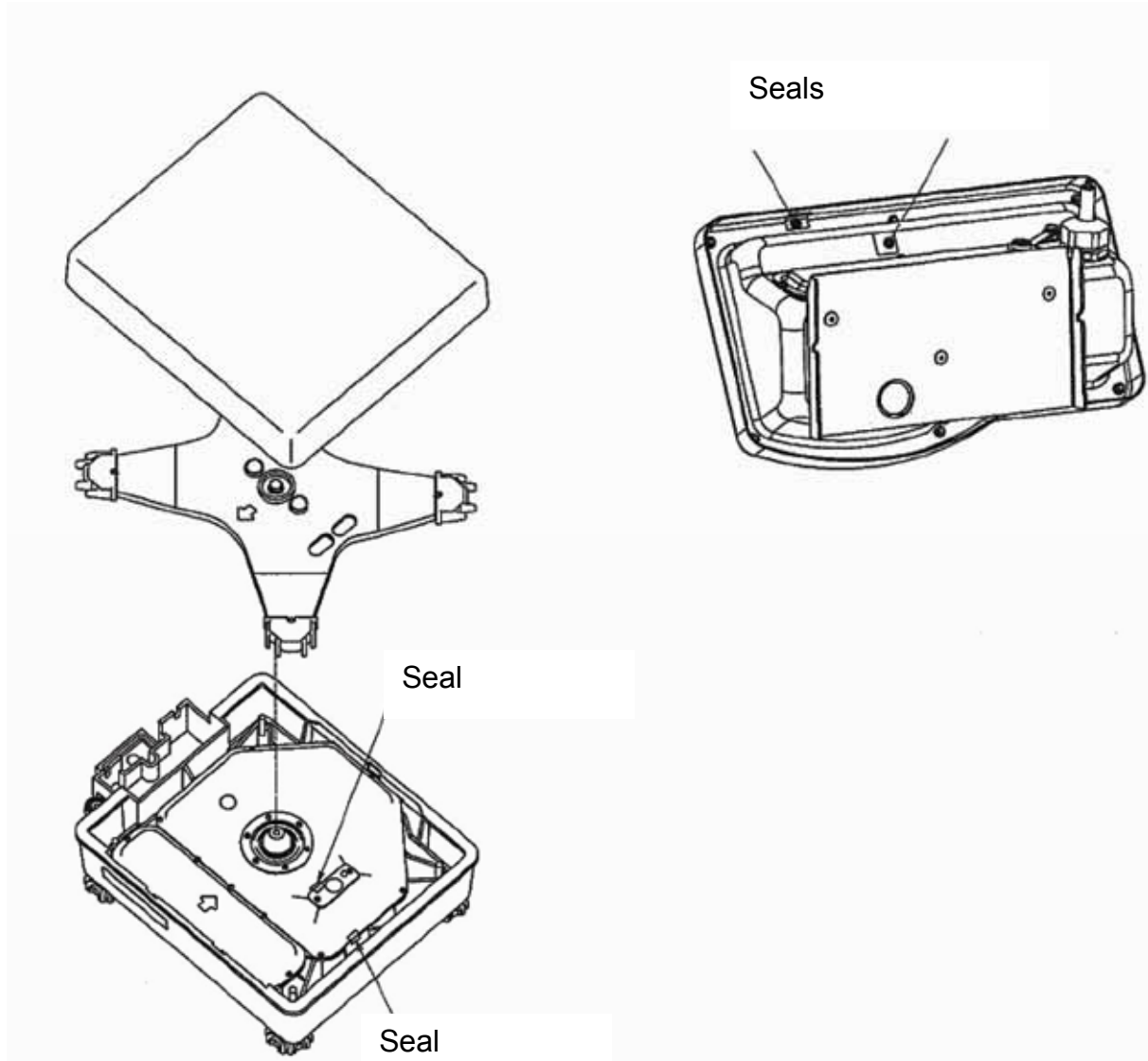
- **Position left: enabled**
- **Position right: Calibration position**

Position calibration switch



After verification the balance is sealed at the indicated positions.
Verification of the balance is invalid without the "seal".

Position of the "official seals":



Balances with obligation to verify must be taken out of operation if:






- **The weighing result of the balance is outside the error limit.** Therefore, in regular intervals load balance with known test weight (ca. 1/3 of the max. load) and compare with displayed value.
- **The reverification deadline has been exceeded.**

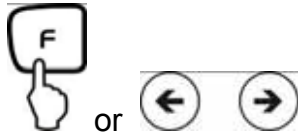
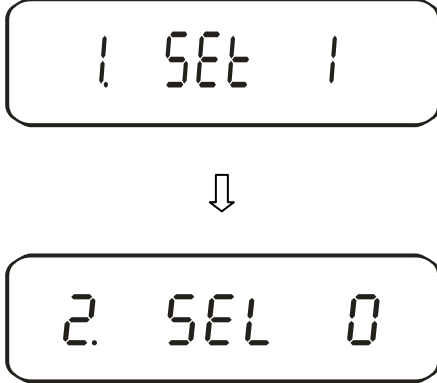
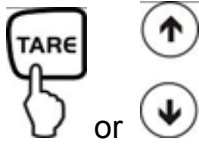
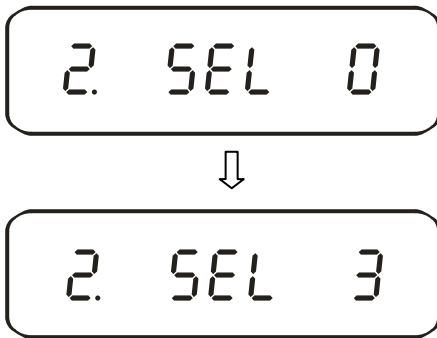


7 Application and configuration menu 1

In the menu the settings of the balance can be modified and functions can be activated. This way, the balance can be adjusted to individual weighing requirements. The menu is structured as follows

- ⇒ **Application menu:** To adjust the balance to user requirements
- ⇒ **Configuration menu 1:** Definition of the basic functions

7.1 User principle of the menu control

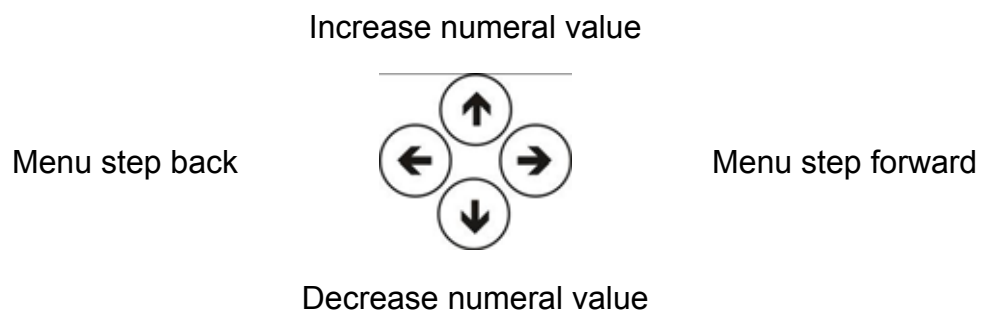
Operation	Display
<p>Switch on balance:</p> 	
<p>Call up menu:</p>  <p>Press for about 4 seconds, until [Func] is displayed.</p>	 <p>When releasing, the first function is displayed [1 SEt 1].</p> 

<p>Changing the function:</p>  <p>Further pressing the keys will call up the various functions of the menu (see table chapter 7.2)</p>	
<p>Changing the parameter:</p>  <p>To change the last digit of the parameter, actuate the TARE-key or the arrow keys.</p>	
<p>Saving your settings:</p>  <p>Leave the function menu and return to weighing mode</p>	

General information about using the arrow keys for entering:







Operation via arrow keys is faster and more comfortable than via the TARE and F key.

Key allocation of the arrow keys:



7.2 Menu overview







The manufacturer's setting has a certain standard configuration. This one is marked with *.

Function	Display  or  	Selectio  or  	Description of the options	
Weighing mode	1 SEt.	* 1	Weighing	
		2	Combination: Weighing/parts counting	
		3	Combination: Weighing/percent determination	
Additional functions	2 SEL	* 0	Off	
		1	Add	
		2	Tolerance weighing	see chapter 12.3, Tab. 1
		3	Combination Tolerance weighing/adding	
Zero balancing	3 RD	0	No zero balancing	
		* 1	Automatic zero balancing is activated.	
Vibration filter	4 Sd.	* 2	Sensitive and fast (very quiet set-up location). ↓	
		3	Robust but slow (very busy set-up location)	
		4		
Display speed	5 rE.	0	Setting for dispensing: Sensitive and fast	
		1		
		2		
		* 3	↓ Insensitive but slow	
Interface (see chpt. 7.2.1)	6 IF.	0	Deactivated	
		* 1	6-digit data format	see chap. 15.4.1
		2	7-digit data format	
		3	extended 7-digit data format	not documented
Adjustment * 1: Factory setting FEJ * 2: Factory setting FES	7 CR.	0	CAL-key turned off	
		* 1	Automatic internal adjustment (only FEJ)	
		2	Adjustment test with internal weight (only FEJ)	
		* 2	External adjustment	
		4	Adjustment test with external weight	

Bar graph	B. bG.	0	Hide bar graph
		* 1	Show bar graph
Automatic turn-off for battery operation (function only exists for battery operation)	9. AP.	0	Automatic turn-off for battery operation (optional) - off.
		* 1	Automatic turn-off for battery operation (optional) - on.
Auto Sleep	A. AS.	0	Off
		* 1	3 minutes after connecting balance to power supply it will go into sleep mode
Units A	b1. uA	* 1	(g)
		2	(kg)
		4	(ct)
Units B	b3. ub	* 0	No unit
		1	(g)
		2	(kg)
		4	(ct)
Display last fractional digit	C. A.1.	0	no
		* 1	Yes; always use this setting!
Multiple range	C. dr	0	no
			Yes
In accordance with ISO/GLP/GMP	E. GLP	* 0	no
		1	Yes
Not documented	E1. out	0	no
		* 1	Yes
	E2. od.	* 0	no
		1	Yes
E3. PF.	* 1	English	
	2	not documented	
Date	F. dAtE	1	Display in year-month-day
		2	Display in month-day-year
		* 3	Display in day-month-year
Time	G. t.o.	* 0	Output - NO
		1	Output - YES
Immediate start	L. dSt.	* 0	When connecting the mains cable, the balance will immediately go into stand-by mode
		1	Balance switches on when plugging in mains power supply
Output interface	n. PrF.	1	not documented
		2	not documented
		* 3	not documented

7.2.1 Parameter for serial interface



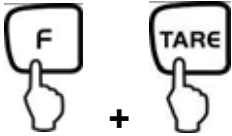


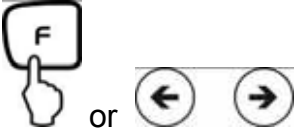


Not shown for menu setting „ $\bar{6}$ 1 F 0“ (interface de-activated).


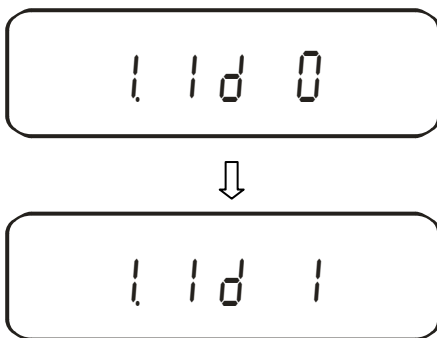


Function	Display  or  	Selectio  or  	Description of the options
Output condition at interface	$\bar{6}$ 1. o.c.	0	No data output
		1	Continuous data output
		2	Continuous data output stable weighing values
		3	Output for stable and instable weighing values after pressing PRINT key
		4	Output for stable weighing value after previous relief of balance
		5	One output for stable weighing value. No output for stable weighing values. Renewed output after stabilization
		6	One output for stable weighing value. Continuous output for instable weighing values.
		* 7	Output of stable weighing values after pressing PRINT key
		8	Single, immediate output after fixed interval (see chpt. 14.5)
		6	Single, immediate output after fixed interval and stable weighing value (see chpt 14.5)
Baud rate	$\bar{6}$ 2. b.L.	* 1	1200 bps
		2	2400 bps
		3	4800 bps
		4	9600 bps
		5	19200 bps

Parity only at setting 6.1F.2 or 6.1F.3	63 PR.	* 0	No parity bit
		1	Odd parity
		2	Even parity
Data Bits only at setting 6.1F.3	64 dL.	* 7	7 bits
		8	8 bits
Stop Bits only at setting 6.1F.3	65 St.	1	1 bit
		* 2	2 bit
not documented	66 uN.	* 0	Always use this setting
		1	
not documented	67 rES.	* 1	Always use this setting
		2	

8 Configuration menu 2



8.1 User principle of the menu control

Operation	Display
<p>Switch on balance:</p> 	
<p>Call up menu:</p>  <p>Hold F-key while TARE key is pressed until [Func 2] is displayed.</p>	 <p>When releasing, the first function is displayed [1. 1d. 0]</p> 
<p>Changing the function:</p>  <p>Further pressing the keys will take you through the various functions of the menu.</p>	 <p style="text-align: center;">↓</p> 

<p>Changing the parameter:</p>  <p>To change the last digit of the parameter actuate TARE-key or arrow keys.</p>	
<p>Saving your settings:</p>  <p>Leave the menu and return to weighing mode</p>	

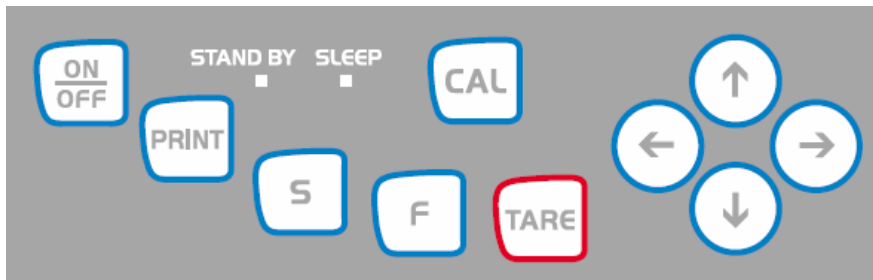
8.2 Menu overview

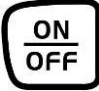





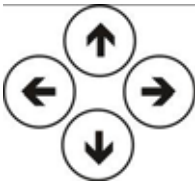


The manufacturer's setting has a certain standard configuration. This one is marked with *.

Function	Display 	Choice 	Description of the options
Setup balance ID no.	1. 1d	*0	Off
		1	On
Overwriting the adjusting weight Caution: Modifications may only be carried out by specialized personnel!	2. rLR	*0	Off
		1	On

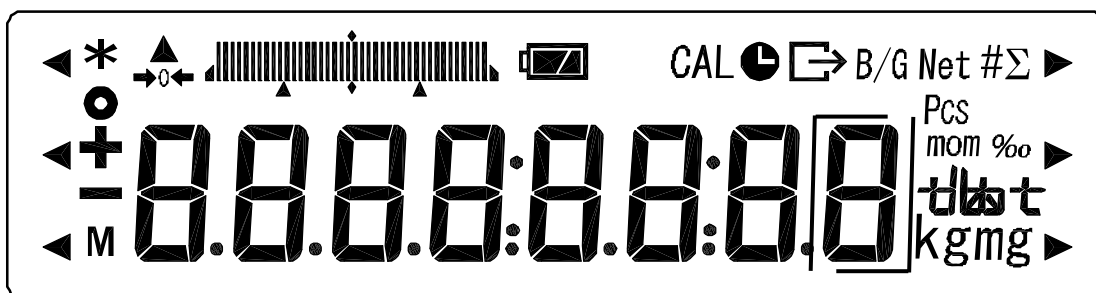
9 Operation

9.1 Keyboard overview



Choice	Function
	<ul style="list-style-type: none"> • Turn on/off
	<ul style="list-style-type: none"> • Output of the weight value on an external device (printer) or PC
	<ul style="list-style-type: none"> • Save function parameters • Addition of displayed values in addition memory • Menu call up "Enter tolerance limits"
	<ul style="list-style-type: none"> • Switching the displayed value (g, ct, Pcs, %) • Entering numeric values • Choosing the function values within the function • Call up individual functions (multiple print) • The entry point will be shifted one spot to the left
	<ul style="list-style-type: none"> • Tare or set weight display to zero • Individual setting within the individual function • Changing the parameters
	<ul style="list-style-type: none"> • Start internal adjustment
	<ul style="list-style-type: none"> • For many entering functions, the arrow keys replace the  or  keys (see chapter 7.1)
LED (green)	<ul style="list-style-type: none"> • "Stand-by" glows if the balance is operated with energy from the power mains but turned off.
LED (red)	<ul style="list-style-type: none"> • "Sleep" has the function of a display saver. It can be deactivated by actuating a key or changing the load.

9.2 Overview of display



Display	Description
g, kg	Gram, Kilogram
→0←	Zeroing display
-	Minus
o	Stability display
Net	Net weight
B/G	Gross /gross weight
Pcs	Parts counting
%	Percent weighing
◀	Tolerance weighing
*	Adding function active
Σ	Total
Ⓛ	Output date/time
M	Balance carries out balance function, e.g. unit count / display of stored value
CAL	Display for adjustment. Signals the adjustment function.
gkg	Weighing unit display
Bar graph	Bar graph
Battery icon	Message for battery mode (optional) see chap. 6.6
□	Display last fractional digit

10 Weighing mode

The manufacturer's standard configuration is "Weighing":

Function [**1 5 E t . 1**], see chapter 7.2.

Under this menu item, you can combine the function "Weighing" with the function "Parts counting" or "Percent determination" (settings see chapter 7.2).




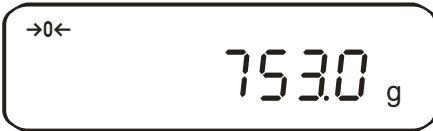

This way, 3 different weighing modes are available for you:

1. Weighing [**1 5 E t . 1**]
2. Weighing/parts counting [**1 5 E t . 2**]
3. Weighing/percent determination [**1 5 E t . 3**]

Apart from selecting a weighing mode, you can activate other functions (Tolerance weighing, adding); see chapter 7. This enables you to display your measuring values according to your requirements.

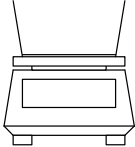





Actuating the **[F]** key will switch the displayed value to the active function (e.g. "g" to "Pcs").

10.1 Weighing

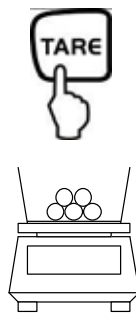

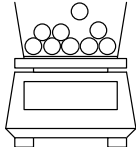
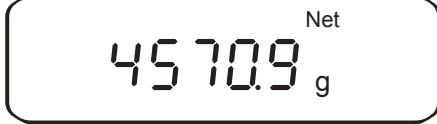
Operation	Display
<p>Switch on balance:</p>  <p>Your balance is ready to weigh as soon as the "0.0" display appears.</p>	<p>The balance will carry out a self-test</p>  <p style="text-align: center;">↓</p> 
<p>Put on items to be weighed, weighed value is displayed.</p>	
 <p>By repeated pressing, switching option of the displayed value into other activated functions/weighing units</p>	

10.1.1 Taring

The dead weight of any weighing container may be tared away by pressing a button, so that the following weighings show the net weight of the goods to be weighed.

Operation	Display
<p>Place empty tare container on the weighing plate. The total weight of the container is displayed.</p> 	
	<p>Reset display to "0":</p>  <p>The weight of the container is now saved internally; in addition the display shows the tare symbol "Net".</p>
<p>Place the goods to be weighed into the tare container.</p> 	<p>Read the weight of the goods on the display.</p> 

The taring process can be repeated any number of times, e.g. when adding several components for a mixture (adding).

 <p>The diagram shows a hand pointing to a button labeled 'TARE' above a balance scale. The scale has a weighing pan containing three small circles representing components.</p>	<p>Reset display to "0":</p>  <p>The digital display shows '0.0' with 'Net' above it and 'g' below it.</p> <p>The total weight of the container is tared away.</p>
<p>Add more components into the weighing container (adding).</p>  <p>The diagram shows the same balance scale as in the first step, but now with five small circles on the weighing pan.</p> <p>Now read off the weight of the added item to be weighed on the display.</p>	 <p>The digital display shows '4570.9' with 'Net' above it and 'g' below it.</p>

NOTE:





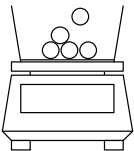



The balance is able to only store one taring value at a time.
When the balance is unloaded the saved taring value is displayed with negative sign.
Remove all items from the weighing plate in order to delete the stored tare value and subsequently press the TARE key.
The taring process can be repeated any number of times. The limit is reached when the whole weighing range is exhausted.


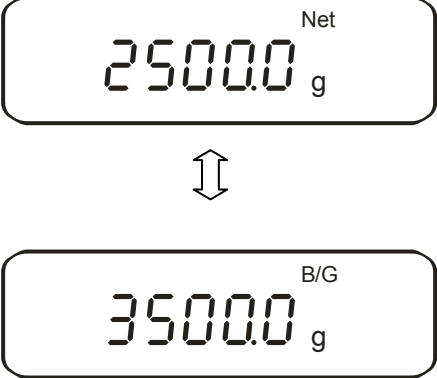
10.1.2 Net/gross

The dead weight of any weighing container may be tared away by pressing a button. For subsequent weighings the net weight of the goods to be weighed as well as the gross weight goods + taring container can be displayed.

Condition:

- Function [! 5E£. !] active (see chapter 7)

Operation	Display
<p>Place empty tare container on the weighing plate. The total weight of the container is displayed.</p> 	
	<p>Reset display to "0":</p>  <p>The weight of the container is internally stored, the display shows the tare symbol "Net".</p>
<p>Place the goods to be weighed into the tare container.</p> 	<p>The net weight of the goods to be weighed is displayed.</p> 
	<p>The gross weight (goods + taring container) is displayed, the display shows the gross symbol "B/G".</p> 

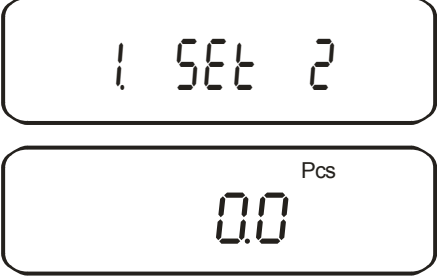

 <p>Use the F key to switch from net weight to gross weight or vice versa</p> <p>This procedure may be repeated any number of times (max. weighing range of the balance).</p>	
--	--



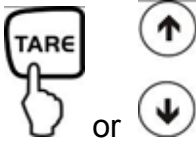
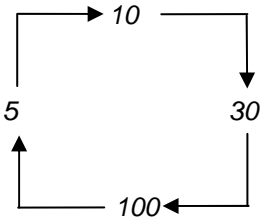
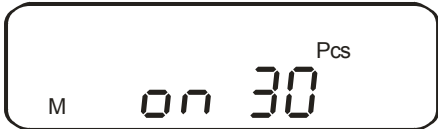



10.2 Weighing/parts counting


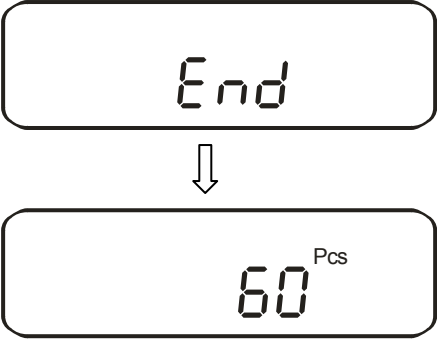
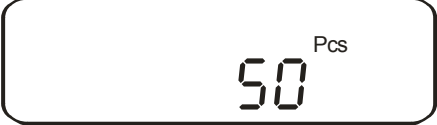

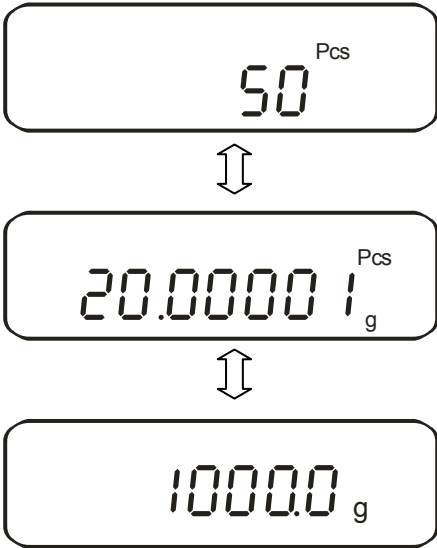
With parts counting you can either count parts into a container or remove parts from a container. To count a greater number of parts the average weight per part has to be determined with a small quantity (reference quantity). The larger the reference quantity, the higher the counting exactness. High reference must be selected for small parts or parts with considerably different sizes.

The process has four steps:

- Tare the weighing container
- Determine the reference unit
- Weigh in the reference unit
- Count the items

Operation	Display
<p>Activate function [1 SET 2] (see chapter 7).</p> <p>The display shows the reference quantity symbol "Pcs".</p>	
 <p>If you are using a weighing container</p>	

<p>Determine the reference unit:</p>  <p>Press for about 4 seconds, until [U. Set.] is displayed, then release</p>	<p>The display shows flashing the last saved reference quantity.</p>  <p>The display e.g. 10 Pcs. prompts you to enter 10 pieces as reference.</p>
<p>Change reference quantity:</p>  <p>Use the TARE-key or the arrow keys to switch between the following reference quantities:</p>  <p>Important: The larger the reference quantity, the more accurate the parts counting.</p>	
<p>Weigh in the reference unit:</p> <p>Place as many parts to count on the balance as the set reference quantity requires.</p> 	<p>The reference quantity is displayed flashing.</p>  <p>The balance offers the possibility of reference optimizing. If this is not desired, press the F key.</p>
<p>Reference optimization:</p> <p>By adding more pieces (up to the 3-fold quantity), you can optimize the reference. At every reference optimization, the reference weight is calculated anew. As the additional pieces increase the base for the calculation, the reference also becomes more exact.</p>	

 <p>The reference weight is saved. Remove reference weight.</p>	
<p>Count the items: Now you can fill the items to be counted into the container. The respective quantity is shown in the display.</p>	
 <p>By repeated pressing, switching option of the displayed value of the displayed value e.g. in:</p> <ul style="list-style-type: none"> • Number of parts placed on balance "Pcs" <ul style="list-style-type: none"> ⇕ • Average part weight "g/Pcs" <ul style="list-style-type: none"> ⇕ • Weight of parts placed on balance in "g" 	

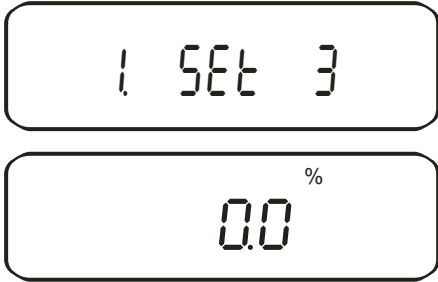



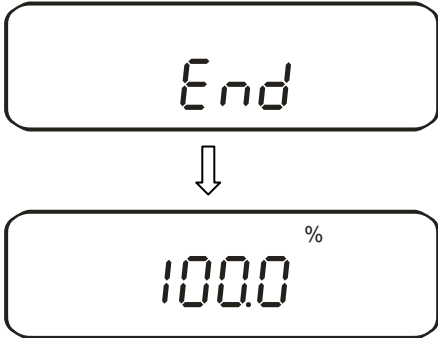
NOTE:

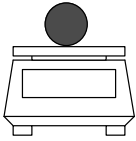



- If the error message "**Sub**" appears, the triple quantity has been exceeded.
- If the error message "**L-Err**" appears the smallest counting weight has not been reached.
- If the "**Add**" error message appears, the applied number of items is too small for correct determination of the reference. For reference, place more parts on the balance.

10.3 Weighing/percent determination

Percent weighing allows to display weight in percent, in relation to a reference weight. The displayed weighing value is stored as a standard percent value (default setting: 100%).

10.3.1 Entering the reference weight by weighing

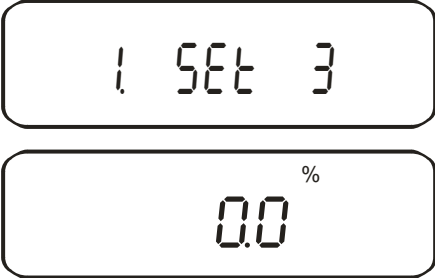

Operation	Display
<p>Activate function [1 5Et 3] (see chap. 7).</p> <p>The display shows the %-symbol.</p>	
<p>Determine reference weight:</p>  <p>Press for about 4 seconds, until [P. 5Et] is displayed, then release</p>	<p>The display shows flashing the last saved reference weight</p>
<p>Put on reference weight (=100 %)</p> 	
 <p>An acoustic signal sounds; the reference weight is saved.</p> <p>Remove reference weight.</p>	



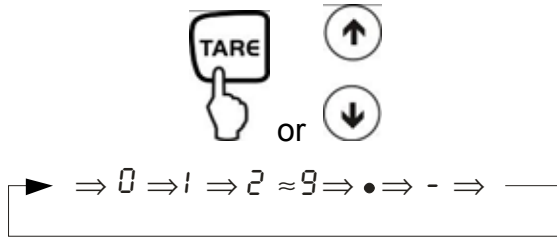
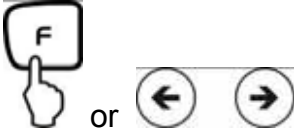

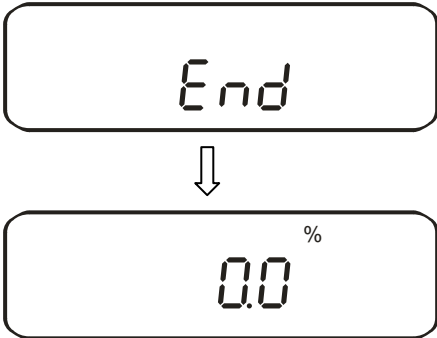
<p>From now, the added weight is shown in %.</p> 	
<p>By repeated pressing, switching option of the displayed value in „g“ or %“</p> 	

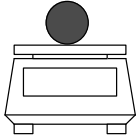



NOTE:

- If the error message “**o-Err**“ is displayed, the reference weight is outside the weighing range
- The 100% reference is preserved until the balance is disconnected from the mains.

10.3.2 Numeric entering of the reference weight

Operation	Display
<p>Activate function [1 5Et 3] (see chap. 7).</p> <p>The display shows the %-symbol.</p>	
<p>Determine reference weight:</p>  <p>Press for about 4 seconds, until [P. 5Et] is displayed, then release</p>	<p>The display shows flashing the last saved reference weight</p>

	 <p>With "0" flashing, you are prompted to enter the reference weight numerically</p>
<p>Entering the numeric value:</p>  <p>Any time you press TARE-key or the arrow key, you will go through the numbers 0-9, decimal dot and minus</p>	
<p>Select the number to be changed (the active position flashes):</p> 	
 <p>An acoustic signal sounds; the entered reference weight is saved</p>	

<p>From now, the added weight is shown in %.</p> 	
<p>By repeated pressing, switching option of the displayed value in „g“ or %“</p> 	

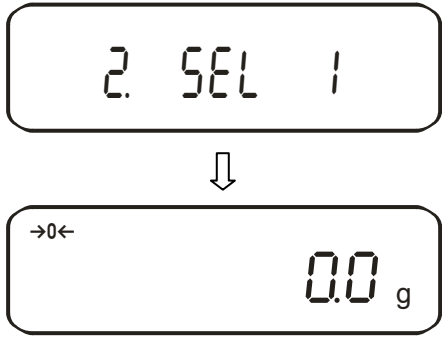




NOTE:


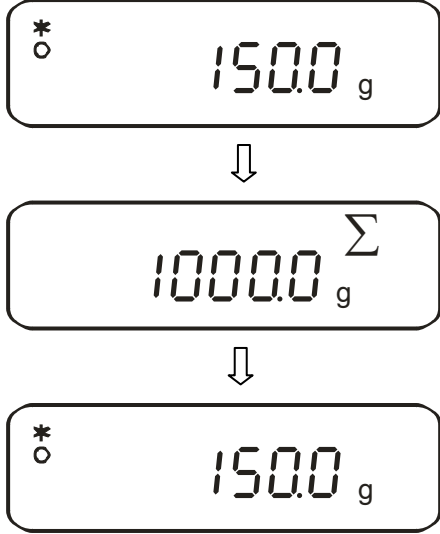


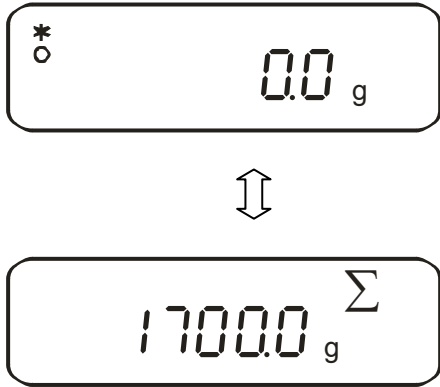


- If the error message “**o-Err**“ is displayed, the reference weight is outside the weighing range
- The 100% reference is preserved until the balance is disconnected from the mains.

11 Adding of displayed values

Any number or individual weighings are automatically added to a total, e.g. all individual weighings of a batch.

The add function is possible in all functions of the weighing mode (Weighing/parts counting/ percent determination).

Operation	Display
1. Activate function [2 SEL 1] (see chap. 7).	
2. Place the weight A , wait until the stability display [O] appears	
3.  The displayed value is added into the total adding memory. The total [Σ] is briefly shown	
4. Remove weight. As soon as the symbol [*] appears further weights can be placed on the balance	
5. Wait until the balance zero display is shown, then place weight B on balance	

<p>6. Wait until stability display is shown [O] :</p>  <p>The displayed value is added into the total adding memory. The total [Σ] is briefly shown</p>	
<p>Remove weight and place further weights on balance; for each weight, repeat step 4 to 6</p>	
<p>7. Total of all individual weighings:</p>  <p>By repeatedly pressing the F-key, switching of the displayed value into other activated functions.</p>	
<p>8. Delete total added memory: Display total (step 7), then press TARE-key.</p> 	

Adding of displayed values is possible without removing the weight. To do so, in step 4 press Tare-key instead of removing the weight.

12 Weighing with tolerance range

12.1 General Information

This balance can be used as dispensing as well as sorting balance; the respective lower tolerance limit as well as upper tolerance limit can be programmed. An acoustic signal supports portioning, dispensing or sorting.

- In the menu, (see chapter 7) activate the tolerance weighing function:

[2.5EL.2]

or the combination tolerance weighing/adding (tolerance control on the respective poured quantity):

[2.5EL.3]

Entering limits is possible for the following functions:

- Weighing
- Parts counting
- Percent determination

There are several functions available for tolerance control of the balance **KERN FEJ**. There are two different ways to carry out evaluation of limits:

1. Evaluation of absolute values **[24. tYP. 1]**:
An exact reference value (e.g. 1 kg) is set.
2. Evaluation with difference values **[24. tYP. 2]**:
An upper limit and a lower limit for a reference value are set.

Example:

	Reference value	Lower limit	Upper limit
Poured quantity	1,000.0 g	970.0 g	1,050.0 g
Evaluation of absolute values	1,000.0 g	970.0 g	1,050.0 g
Evaluation with difference values	1,000.0 g	-30.0 g	50.0 g

There are two different ways to set the tolerance limits:

1. Place the values (object) on the balance -
 > Save this value
2. Numeric entering of values -
 > Enter the limits via keyboard.

NOTE:

- ⇒ If a limiting value was set it remains saved until the balance is turned off.
- ⇒ For the functions weighing, counting, percent individual limits can be set.
- ⇒ When entering the limits please pay attention to the type of evaluation that was set.

12.2 Display of the results

12.2.1 For 2 limits

The triangular tolerance marker (◀) in the upper part of the display shows whether the goods to be weighed are within the two tolerance limits. The tolerance marker is only in operation during operating mode tolerance weighing; it is otherwise not visible.

The tolerance marker provides the following information:



Goods to be weighed above tolerance limit



Goods to be weighed within tolerance range

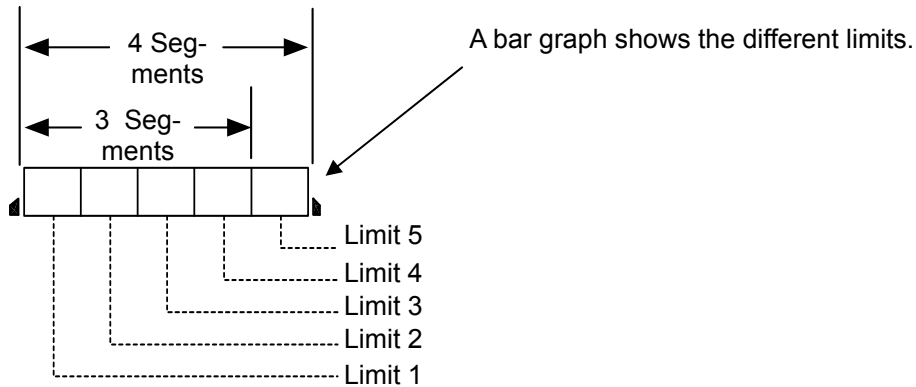


Goods to be weighed below tolerance limit











Display Result	If a point is set as lower limit	If two points are set as upper and lower limit
+ (high)	No display	Weight > Upper limit
TOL ✓ (OK)	Lower limit ≤ Weight	Lower limit ≤ Weight ≤ Upper limit
- (low)	Lower limit > Weight	Lower limit > Weight

12.2.2 For 3 or 4 limits







If 3 or 4 limiting points are to be set, this is displayed in the bar graph. The length of the represented bars indicates where within the tolerance range the weighed of the goods to be weighed are.



12.3 Basic settings for weighings with tolerance range

Operation	Display
1. Activate tolerance weighing function [2.5EL.2] or [2.5EL.3] (see chap. 7).	
2. Selection of tolerance parameters  or   <p>Any time you press the F-key you can select between the following settings, see tab. 1.</p>	 <p>The first parameter for setting the tolerance marker appears.</p>
3. Changing the parameter value  or  	 <p style="text-align: center;">↓</p> 

Tab. 1:








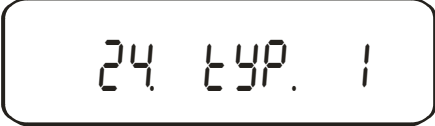


Function	Display  or  	Selectio  or  	Description of options
Display conditions of the tolerance marker	21. Co.	* 1	Tolerance marker is always displayed, even if standstill control is not yet displayed.
		2	Tolerance marker is only displayed in connection with standstill control.
Tolerance range	22. Li.	0	Tolerance marker is only displayed above zero range (mind + 5).
		* 1	Tolerance marker is displayed for the whole range.
Number of limiting points	23. Pi	1	1- Limiting point (OK/ -)
		*2	2- Limiting points (+/OK/-)
		3	3- Limiting points (1-4)
		4	4- Limiting points (1-5)
Assessment	24. tYP	* 1	Evaluation for absolute values
		2	Evaluation for difference values (with reference weight)
Signal at limit 1	25. bu1	*0	No signal at limit 1(-)
		1	Signal at limit 1 (-)
Signal at limit 2	26. bu2	*0	No signal at limit 2(Ok)
		1	Signal at limit 2(Ok)
Signal at limit 3	27. bu3	*0	No signal at limit 3(+)
		1	Signal at limit 3(+)
Signal at limit 4	28. bu4	*0	No signal at limit
		1	Signal at limit 4
Signal at limit 5	29. bu5	*0	No signal at limit 5
		1	Signal at limit 5
Display of Results	2A. LG	* 1	Display via +, OK or -
		2	For setting 2 limits display in bar graph is possible
Output setting	2b r.o.c.		Not documented


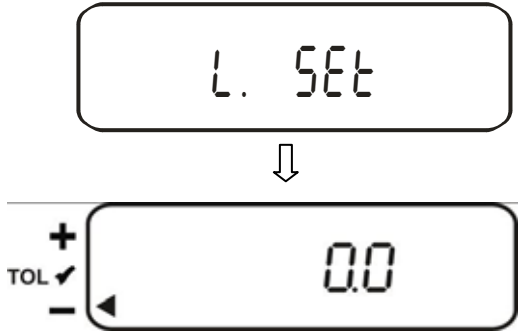
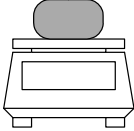

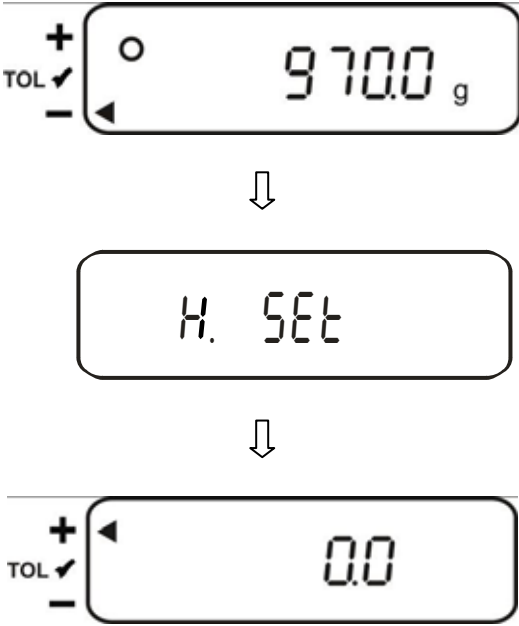
12.4 Evaluation of absolute values

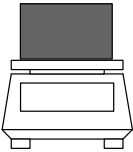

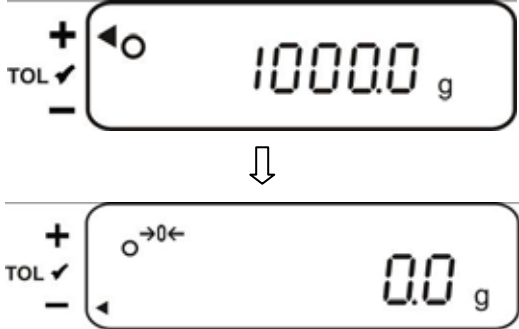
12.4.1 Entering 2 limits by weighing

Important information!

Always begin by entering the lower limit value, followed by the upper limit value. Enter.


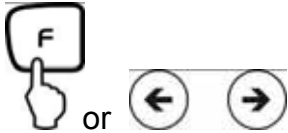


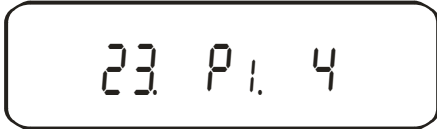
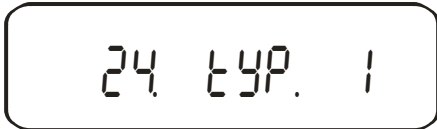

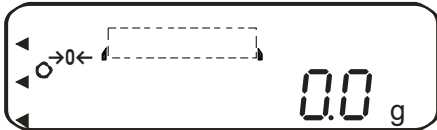
Operation	Display
1. Activate tolerance weighing function [2.5EL.2] or [2.5EL.3] (see chap. 7).	 
2. Actuate required parameter selection  or  until [23. P1.1] or [24. tYP.1] is displayed; more settings of your choice (see tab.1, chapter 12.3) are carried out in the same way	  Parameter selection for 2 limiting points:  Parameter selection for absolute value: 
3. Leave function menu 	 The balance is now in tolerance weighing mode; the tolerance marker (◀) appears

<p>4. Entering limiting values:</p>  <p>Press for about 4 seconds, until [L. SEt] is displayed, then release</p>	 <p>The flashing display (last saved value) prompts you to enter the lower limiting value (L. SEt)</p>
<p>5. Place sample for the lower (i.e. smaller) limiting value on the weighing plate:</p> 	
<p>6. Save:</p> 	<p>An acoustic signal sounds, the saved lower limit is briefly displayed.*</p>  <p>The flashing display (last saved value) prompts you to enter the upper limiting value (H. SEt)</p>

<p>7. Place sample for the upper (i.e. larger) limiting value on the weighing plate:</p> 	
<p>8. Save:</p>  <p>The balance returns to tolerance weighing mode. From here evaluation takes place whether the goods to be weighed are within the two tolerance limits.</p>	<p>An acoustic signal sounds, the saved upper limit is briefly displayed.</p> 

* If you want to set for your tolerance weighing only one limit point (parameter selection [23. P 1. 1]), ignore step 7 and 8.

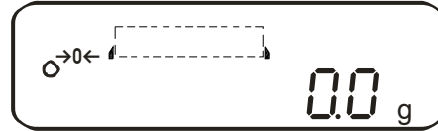
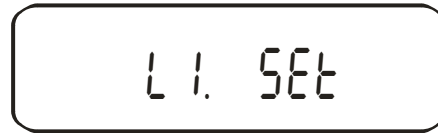
12.4.2 Entering 3 or 4 limits by weighing

Operation	Display
<p>1. Activate tolerance weighing function [2.5EL.2] or [2.5EL.3] (see chap. 7).</p>	 <p style="text-align: center;">↓</p>
<p>2. Actuate required parameter selection</p>  <p style="text-align: center;">or</p> <p style="text-align: center;">until</p> <p>[23. Pl. 1] or [24. tYP. 1] is displayed;</p> <p>more settings of your choice (see chap. 12.3) are carried out in the same way</p>	 <p style="text-align: center;">↓</p> <p>Parameter selection for 3 limiting points:</p>  <p>Parameter selection for 4 limiting points:</p>  <p>Parameter selection for absolute value:</p> 
<p>3. Leave function menu</p> 	 <p>The balance is now in tolerance weighing mode; as a tolerance marker, the bar graph is displayed.</p>

4. Entering limiting values:

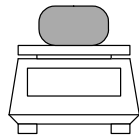


Press for about 4 seconds, until [L 1 SEt] is displayed, then release



The flashing display (last saved value) prompts you to enter the first lower limiting value (L1 SEt)

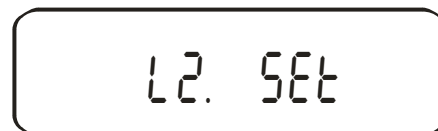
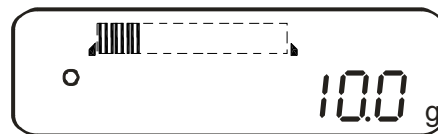
5. Place sample for the first limiting value on the weighing plate:



6. Save:

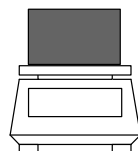



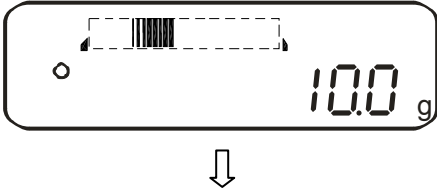
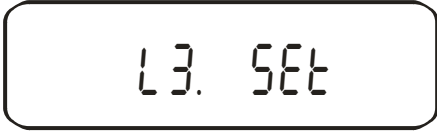

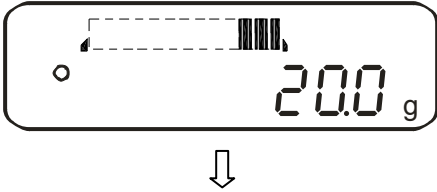
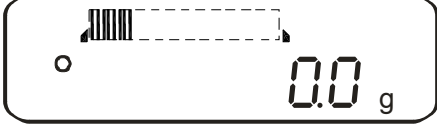
An acoustic signal sounds, the saved first limiting value is briefly displayed.*






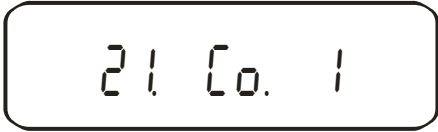
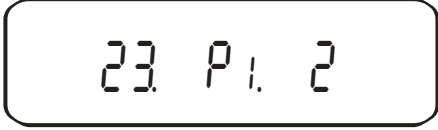




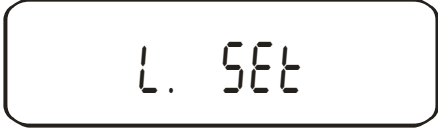
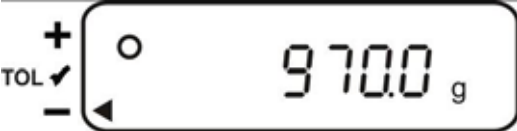
The flashing display (last saved value) prompts you to enter the second limiting value (L2 SEt)



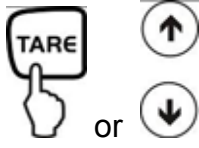
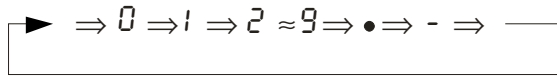
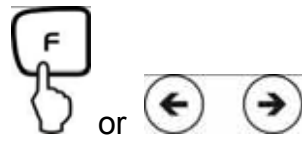
7. Place sample for the second limiting value on the weighing plate:


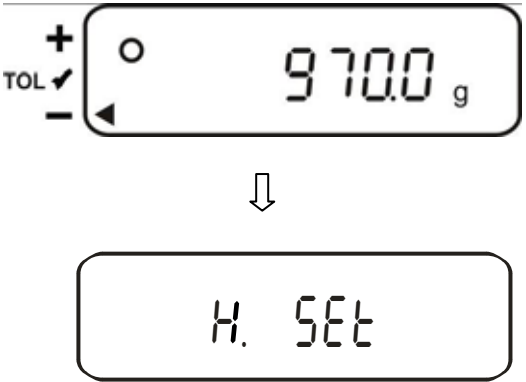

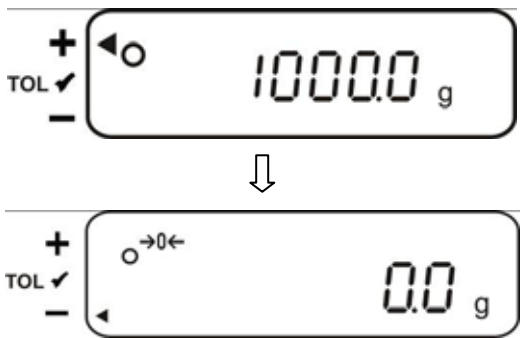


<p>8. Save:</p> 	<p>An acoustic signal sounds, the saved second weighing value is briefly displayed.</p>   <p>The flashing display (last saved value) prompts you to enter the third limiting value (L 3.5Et)</p>
<p>9. To enter 3rd and 4th limiting value, repeat steps 7 and 8</p>	
<p>10. Save:</p>  <p>The balance returns to tolerance weighing mode. From here, evaluation is carried out whether the goods to weighed are within the tolerance limits.</p>	<p>An acoustic signal sounds, the last saved 3rd or 4th limiting value is briefly displayed.</p>  

12.4.3 Numeric entering of 2 limits

Operation	Display
<p>1. Activate tolerance weighing function [2.5EL.2] or [2.5EL.3] (see chap. 7).</p>	 <p style="text-align: center;">↓</p>
<p>2. Actuate required parameter selection</p>  <p style="text-align: center;">or</p>  <p style="text-align: center;">until</p> <p>[23. Pl. 1] or [24. tYP. 1] is displayed;</p> <p>more settings of your choice (see chap. 12.3) are carried out in the same way</p>	 <p style="text-align: center;">↓</p> <p>Parameter selection for 2 limiting points:</p>  <p>Parameter selection for absolute value:</p> 
<p>3. Leave function menu</p> 	 <p>The balance is now in tolerance weighing mode; the tolerance marker (◀) appears</p>
<p>4. Entering limiting values:</p>  <p>Press for about 4 seconds, until [L. 5EL] is displayed, then release</p>	 <p style="text-align: center;">↓</p>  <p>The last saved limiting value is displayed flashing</p>

<p>5.</p> 	<p>Display changes to flashing "zero"</p>  <p>The flashing display prompts you to numerically enter the lower limit</p>
<p>6. Entering numeric value for the lower limit</p>   <p>Any time you press the TARE-key or the arrow keys you will go through the numbers 0-9, decimal dot and minus</p>	
<p>Select the number to be changed (the respective active position flashes)</p> 	

<p>7. Save:</p> 	<p>An acoustic signal sounds, the saved lower weighing value is briefly displayed.</p>  <p>The flashing display (last saved value) prompts you to enter the upper limiting value</p>
<p>8. To enter the numeric value for the upper limiting value, repeat steps 5 - 6</p>	
<p>9. Save:</p>  <p>The balance returns to tolerance weighing mode. From here evaluation takes place whether the goods to be weighed are within the two tolerance limits.</p>	<p>An acoustic signal sounds, the saved upper limit is briefly displayed.</p> 





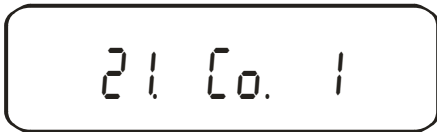

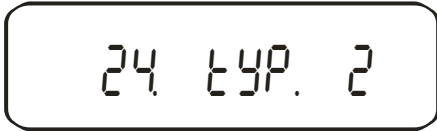


To enter 3 or 4 limiting values [L 1 SEt] - [L 3 SEt] or [L 4 SEt], repeat steps 5 to 7 (see also chapter 12.4.2).


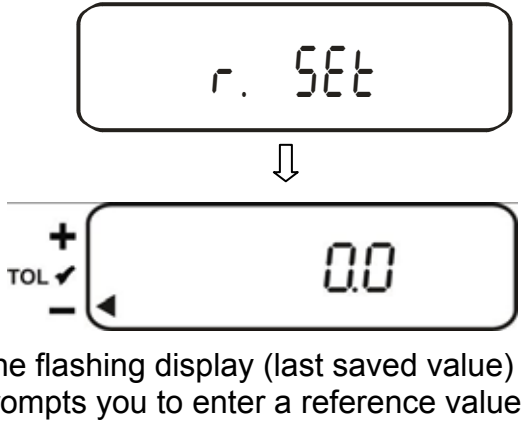
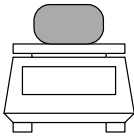

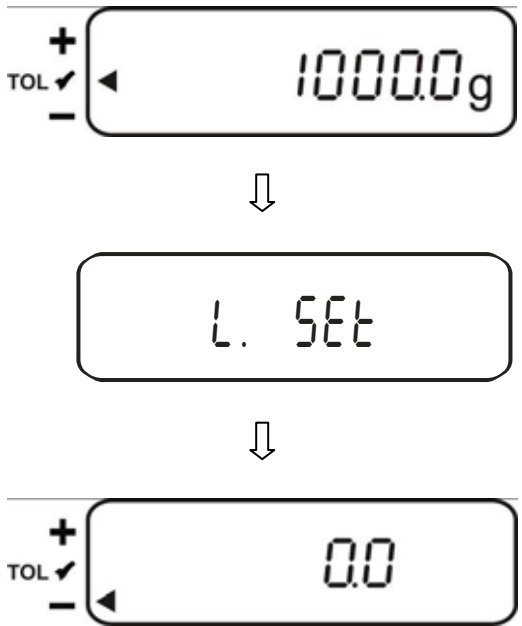
12.5 Evaluation with difference values

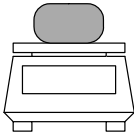


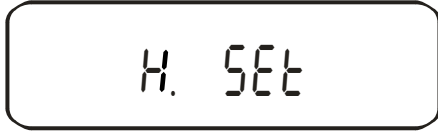
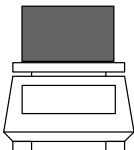

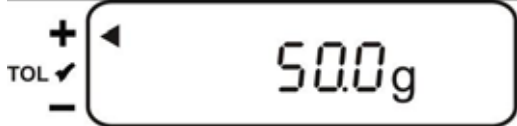

12.5.1 Entering 2 limits by weighing

Important information!

Always begin by entering the lower limit value, followed by the upper limit value. Enter.

Operation	Display
<p>1. Tolerance weighing function [2.5EL.2] or [2.5EL.3] (see chapter 7).</p>	 <p style="text-align: center;">↓</p>
<p>2. Actuate required parameter selection</p>  <p style="text-align: center;">or</p>   <p style="text-align: center;">until</p> <p>[23. P1.2] or [24. tYP.2] appears; more settings of your choice (see chap. 12.3) are carried out in the same way</p>	 <p style="text-align: center;">↓</p> <p>Parameter selection for 2 limiting points:</p>  <p>Parameter selection for difference value:</p> 
<p>3. Leave function menu</p> 	 <p>The balance is now in tolerance weighing mode; the tolerance marker (◀) appears</p>

<p>4. Entering a reference value:</p>  <p>Press for about 4 seconds, until [r.5Et] appears, then release.</p>	 <p>The flashing display (last saved value) prompts you to enter a reference value</p>
<p>5. Place reference weight onto weighing plate:</p> 	
<p>6. Save</p> 	<p>An acoustic signal sounds, the saved reference value is briefly displayed.*</p>  <p>The flashing display (last saved value) prompts you to enter the lower limiting value</p>


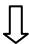


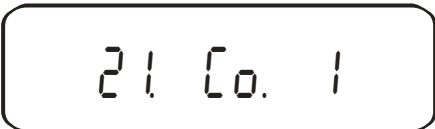


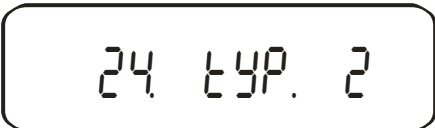


<p>7. Place sample for the first limiting value on the weighing plate:</p> 	
<p>8. Save</p> 	<p>An acoustic signal sounds, the saved lower difference value is briefly displayed.</p>  <p style="text-align: center;">↓</p>  <p>The flashing display (last saved value) prompts you to enter the upper limiting value</p>
<p>9. Place sample for the upper (i.e. larger) limiting value on the weighing plate:</p> 	
<p>10. Save</p>  <p>Remove sample from weighing plate. The balance returns to tolerance weighing mode. From here evaluation takes place whether the goods to be weighed are within the two tolerance limits.</p>	<p>An acoustic signal sounds, the saved upper difference value is briefly displayed.</p>  <p style="text-align: center;">↓</p> 


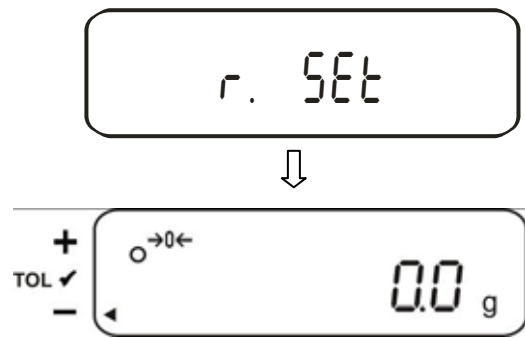


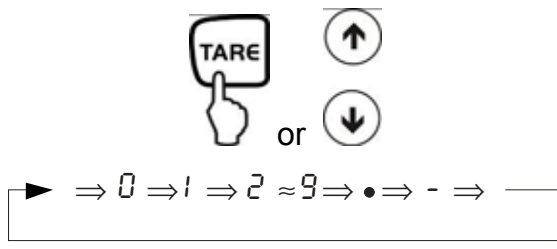
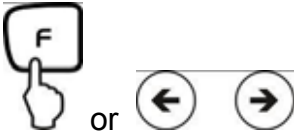
* If you want to set for your tolerance weighing only one limit point (parameter selection [23. P i. 1]), the input herewith is finished.


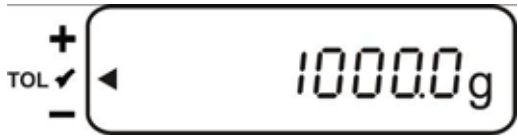
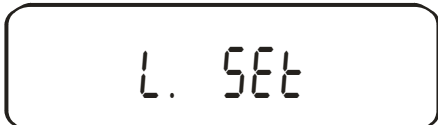



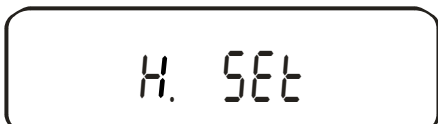
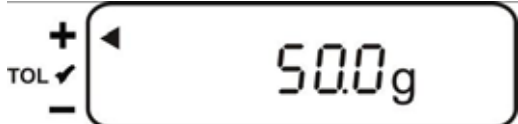
12.5.2 Entering 3 or 4 limits by weighing


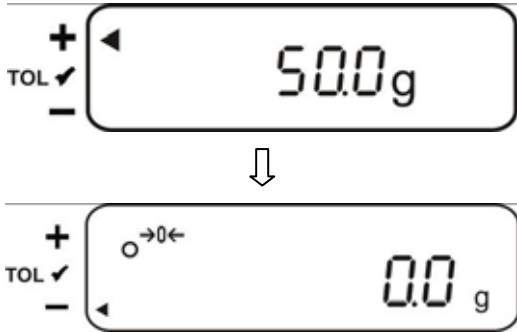
To enter 3 or 4 limiting values [L 1 SEt] - [L 3 SEt] or [L 4 SEt], repeat steps 7 and 8 (see also chapter 12.4.2).

12.5.3 Numeric entering of 2 limits

Operation	Display
1. Activate tolerance weighing function [2.5EL.2] or [2.5EL.3] (see chap. 7).	 
2. Actuate required parameter selection  or  until [23. P1.2] or [24. tYP.2] appears; more settings of your choice (see chap. 12.3) are carried out in the same way	  Parameter selection for 2 limiting points:  Parameter selection for difference value: 
3. Leave function menu 	 The balance is now in tolerance weighing mode; the tolerance marker (◀) appears


<p>4. Entering a reference value:</p>  <p>Press for about 4 seconds, until [r.5Et] appears, then release.</p>	 <p>The last saved reference weight appears flashing</p>
<p>5.</p> 	<p>Display changes to flashing "zero"</p>  <p>The flashing display prompts you to numerically enter a reference weight</p>
<p>6. Entering a numeric value</p>  <p>Any time you press TARE-key or the arrow key, you will go through the numbers 0-9, decimal dot and minus</p>	
<p>Select the number to be changed (the respective active position flashes)</p> 	

<p>7. Confirm</p> 	<p>An acoustic signal sounds, the saved reference weight is briefly displayed.</p>  <p style="text-align: center;">↓</p>  <p>The flashing display (last saved value) prompts you to enter the lower difference value</p>
<p>8. Entering lower limit Repeat steps 5 and 6</p>	
<p>9. Confirm</p> 	<p>An acoustic signal sounds, the saved lower difference value is briefly displayed.</p>  <p style="text-align: center;">↓</p>  <p>The flashing display (last saved value) prompts you to enter the upper difference value</p>
<p>10. Entering upper limit Repeat steps 5 and 6</p>	


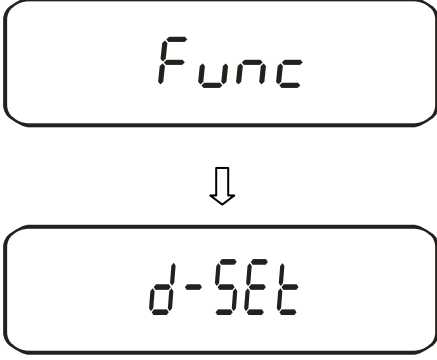
<p>11. Save</p>  <p>The balance returns to tolerance weighing mode. From here evaluation takes place whether the goods to be weighed are within the two tolerance limits.</p>	<p>An acoustic signal sounds, the saved upper difference value is briefly displayed.</p> 
--	---








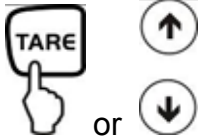
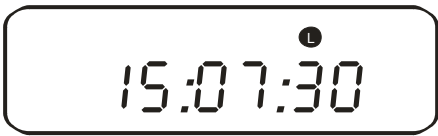




To enter 3 or 4 limiting values [L 1 SEt] - [L 3 SEt] or [L 4 SEt], repeat steps 8 and 9 (see also chapter 12.4.2).

13 Setting date and time

Display symbol []

13.1 Time



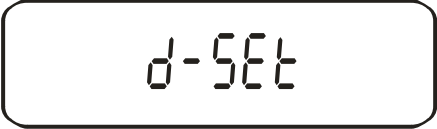




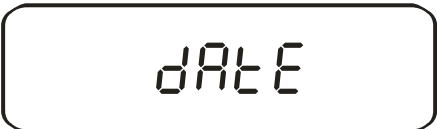

Operation	Display
<p>1. Call up menu</p>  <p>Keep pressed until [d-SEt] appears.</p>	


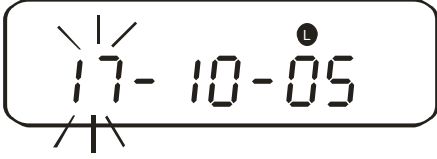

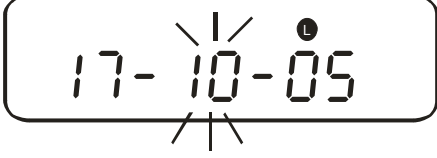
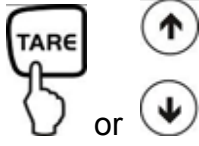



 <p>Press again</p>	 <p style="text-align: center;">↓</p>  <p>Last stored time appears.*</p>
<p>2. Resetting time</p> 	 <p>Time to be changed is flashing</p>
<p>Select the number to be changed (the active position flashes):</p> 	
<p>Changing the numerical value</p> 	
<p>3. Save</p> 	<p>After storing your setting, the date display will appear</p> 
<p>4. Return to weighing mode</p> 	

*Notice: The display can be rounded up (from 30 s) or down (to 29 s) by pressing the TARE-key.

13.2 Date

You can set the display of your data output under menu item [*F. dAtE*] (see menu overview chap. 7.2.).

Operation	Display
<p>1. Call up menu</p>  <p>Keep pressed until [<i>d-SEt</i>] appears.</p>	 <p style="text-align: center;">↓</p> 
 <p>Press again</p>	 <p style="text-align: center;">↓</p>  <p>Last stored time appears</p>
 <p>Press again</p>	 <p style="text-align: center;">↓</p>  <p>Last stored date appears</p>

<p>2. Resetting the date</p> 	 <p>Time to be changed is flashing</p>
<p>Select the number to be changed (the active position flashes):</p> 	
<p>Changing the numerical value</p> 	
<p>3. Save</p> 	<p>After storing your settings, the balance will automatically return to weighing mode.</p> 

14 Description of individual functions

14.1 Auto sleep function

In this function the display of the balance is turned off after 3 minutes without load change or operation to save energy.

Parameter selection: [*A*. *AS* *!*]

The Auto sleep function is indicated by a red LED.

NOTE:

This function is only available when connected to the mains power.

14.2 Auto OFF function

In this function the balance is turned off after 3 minutes without load change or operation to save the battery.


Parameter selection: [*9*. *AP* *!*]

The Auto OFF function is indicated by a green LED.

NOTE:

This function is only available for battery operation.

14.3 Set units function

With this setting you can set different display units (A or B) for one weighing value. Press the -key to choose between units A and B.

14.4 Multiple range display


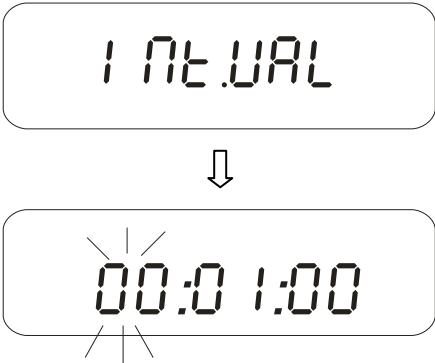


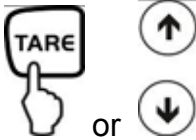



The models FES/FEJ62K0.1DIPM are by standard set as multiple range balance 6.2kg = 0.1g and 62kg= 1g readability. The balance can also be set as single range balance with a readability of 1g.

Parameter selection: [*L*. *dr* *!*]







14.5 Interval output function

This menu item allows you to determine after which interval you wish data output to be carried out. To achieve this, activate the [*5 I. o. c A*] or [*5 I. o. c B*] function in the menu (see chpt. 7.2.1)

14.5.1 Interval setting

Operation	Display
<p>1. Call up menu</p>  <p>Keep pressed until [<i>5 I. o. c A</i>] appears.</p>	 <p>Time to be changed is flashing</p>
<p>2. Set interval</p> <p>Select the number to be changed (the active position flashes):</p> 	
<p>Changing the numerical value</p> 	
<p>3. Save:</p> 	<p>After storing your settings, the balance will automatically return to weighing mode.</p> 

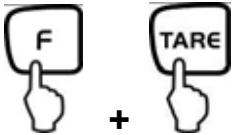


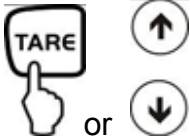





14.5.2 Start/Stop interval output

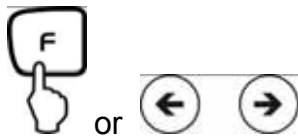

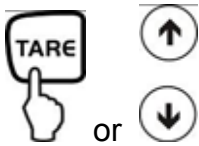





Operation	Display
 <p>Start output</p>	 <p style="text-align: center;">↓</p> 
 <p>Stop output</p>	 <p style="text-align: center;">↓</p>  <p>The balance returns automatically into weighing mode.</p>

14.6 Input balance ID-no.

Display symbol [◀] and [▲]

You can enter a 6-digit number using the characters [0-9], [A-F] and [-]. Space character is displayed as [_].

Operation	Display
<p>1. Call up menu</p>  <p>Hold F-key while TARE –key is pressed until [Func 2] is displayed. see chapter 8.</p>	 <p>When releasing, the first function is displayed [1 . 1 d. 0]</p> 
<p>2. Activate function</p> 	
<p>3. Display serial number</p> 	 <p>The last stored number appears</p>
<p>4. Input serial number</p> 	 <p>Time to be changed is flashing</p>

<p>Select the number to be changed (the active position flashes):</p> 	
<p>Changing the numerical value</p> 	
<p>5. Save:</p> 	<p>Your setting will be stored and the next menu item will appear.</p> 
<p>6. Return to weighing mode</p> 	

15 Data output

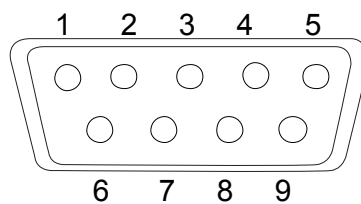
The regular equipment of the balance includes an RS 232C interface and a printer interface.

15.1 RS 232C interface

The RS 232C interface allows a bi-directional data exchange from the FEJ to external devices. This data exchange is asynchronous using ASCII - Code.

Pin allocation of balance output plug:

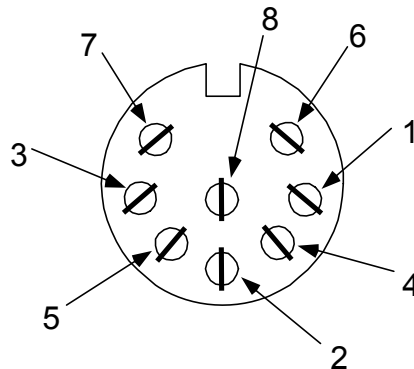
Pin nr.	Signal	Input/Output	Function
1	-		
2	RXD	Input	Receive data
3	TXD	Output	Transmit data
4	DTR	Output	HIGH
5	GND	-	Signal ground
6	-	-	
7	-	-	
8	-	-	
9	-	-	



15.2 Printer interface (unidirectional data exchange)

Pin allocation of balance output plug:

Pin nr.	Signal	Input/Output	Function
1	EXT.TARE	Input	External tare function
2	-		
3	-		
4	TXD	Output	Transmit data
5	GND	-	Signal ground
6	-	-	
7	-	-	
8	-	-	



15.3 Description of interface

The selection of a certain operating mode allows you to set the output format, the output control, the transmission speed and the parity bit. The different options are described in **chap. 7.2.1** under "parameter for serial interface".

15.4 Data output

15.4.1 Format for data transmission

Any of the data formats below may be set by selecting the relevant function on the balance (see menu overview chap. 7.2):

- **6-digit data format**

Consisting of 14 words, including final character; CR=0DH, LF=0AH (CR=balance reverse motion / LF=line feed)

1	2	3	4	5	6	7	8	9	10	11	12	13	14
P1	D1	D2	D3	D4	D5	D6	D7	U1	U2	S1	S2	CR	LF

- **7-digit data format**

1	2	3	4	5	6	7	8	9	10	11	12	13	14	15
P1	D1	D2	D3	D4	D5	D6	D7	D8	U1	U2	S1	S2	CR	LF

NOTE: The 7-digit format is identical to the 6-digit, with the exception of the additional character D8.

- **Extended 7-digit data format**

Not documented

For models e (calibration) = 10 d (legibility) the formats shown below are available:

- **6-digit data format**

1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	
P1	D1	D2	D3	D4	D5	D6	D7	D8	U1	U2	S1	S2	CR	LF	

Consisting of 15 words, including final character; CR=0DH, LF=0AH (CR=balance reverse motion / LF=line feed). Slash "/" is inserted after the e-value.

- **7-digit data format**

1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16
P1	D1	D2	D3	D4	D5	D6	D7	D8	D9	U1	U2	S1	S2	CR	LF

Hinweis: The 7-digit format is identical to the 6-digit format with the exception of the additional character D9.

15.4.2 Signs

P 1 = 1 word

P1	Code	Meaning
+	2 B H	Data is 0 or positive
-	2 D H	Data is negative

15.4.3 Data

6-digit data format (D1-D7): 7 words
 7-digit data format (D1-D8): 8 words
 6-digit data format e = 10 d (D1-D7): 7 words
 7-digit data format e = 10 d (D1-D8): 8 words

D1-D7, D8, D9	Code	Meaning
0 - 9	30 H – 39 H	Data 0 to 9 (max. 6 characters in 6-digit format)
•	2 EH	Decimal point, position not fixed
Sp	20 H	Space character, leading zero suppressed
/	2 FH	Slash "/" is inserted after the e-value.

15.4.4 Units

U 1, U 2 = 2 words as ASCII-Codes

U1	U2	Code		Meaning	Symbol
(SP)	G	20H	47H	Gram	g
K	G	4BH	47H	Kilogram	kg
C	T	43H	54H	Karat	ct
P	C	50H	43H	Quantity	Pcs
(SP)	%	50H	25H	Percent	%

15.4.5 Result evaluation for balances with tolerance range

S 1 = 1 word

S1	Code	Meaning	
L	4CH	Goods to be weighed below tolerance limit	1- or 2 end points
G	47H	Goods to be weighed within tolerance range	
H	48H	Goods to be weighed above tolerance limit	
1	31H	Limit 1	3- or 4 end points
2	32H	Limit 2	
3	33H	Limit 3	
4	34H	Limit 4	
5	35H	Limit 5	
T	54H	Value total	File type
U	55H	Weight value	
(SP)	20H	No rating	
d	64H	Gross	

15.4.6 Data status

S 2 = 1 word

S 2	Code	Meaning
S	53 H	Data stabilized *
U	55 H	Data not stabilized (fluctuating) *
E	45 H	Data error, all data apart from S 2 unreliable Balance indicating error (o-Err, u-Err)
sp	20 H	No special status

15.4.7 Interval data output

When an interval output is started or stopped, this will result in the output of a header or footer line.

Header

- consisting of 15 words

1	2	3	4	5	6	7	8	9	10	11	12	13	14	15
-	-	-	-	-	-	-	-	-	-	-	-	-	-	-

Footers

- Two line feeds are inserted.

15.4.8 Output time

1	2	3	4	5	6	7	8
h	h	:	m	m	:	s	s

* hh: Hours (00-23), mm: Minutes (00-59) min: Seconds (00-59) s:

15.5 Remote control instructions

C1	C2	Code		Meaning
0	0	4FH	30H	No data output
0	1	4FH	31H	Continuous data output
0	2	4FH	32H	Continuous data output stable weighing values
0	3	4FH	33H	Output for stable and instable weighing values after pressing PRINT key
0	4	4FH	34H	Output for stable weighing value after previous relief of balance
0	5	4FH	35H	One output for stable weighing value. No output for stable weighing values. Renewed output after stabilization
0	6	4FH	36H	One output for stable weighing value. Continuous output for instable weighing values.
0	7	4FH	37H	Output of stable weighing values after pressing PRINT key
0	8	4FH	38H	Single immediate output
0	9	4FH	39H	Single output after stabilization
0	A	4FH	41H	Single, immediate output after fixed interval (see chap. 14.5)
0	B	4FH	42H	Single, immediate output after fixed interval and stable weighing value (see chap 14.5)

16 Service, maintenance, disposal

16.1 Cleaning

Before cleaning, please disconnect the appliance from the operating voltage.

Please do not use aggressive cleaning agents (solvents or similar agents), but a cloth dampened with mild soap suds.

The weighing terminal has a **pressure compensation device**.

This is underneath the terminal and made of a glued on membrane.

When cleaning please ensure that the **membrane is not damaged** or soiled.

16.2 Service, maintenance

The appliance may only be opened by trained service technicians who are authorized by KERN.

Before opening, disconnect from power supply.

16.3 Disposal

Disposal of packaging and appliance must be carried out by operator according to valid national or regional law of the location where the appliance is used.

17 Instant help

In case of an error in the program process, briefly turn off the balance and disconnect from power supply. The weighing process must then be restarted from the beginning.

Help:

Fault	Possible cause
The displayed weight does not glow.	<ul style="list-style-type: none">• The balance is not switched on.• The mains supply connection has been interrupted (mains cable not plugged in/faulty).• Power supply interrupted.
The displayed weight is permanently changing	<ul style="list-style-type: none">• Draught/air movement• Table/floor vibrations• The weighing plate is in contact with foreign matter.• Electromagnetic fields / static charging (choose different location/switch off interfering device if possible)
The weighing value is obviously wrong	<ul style="list-style-type: none">• The display of the balance is not at zero• Adjustment is no longer correct.• Great fluctuations in temperature.• Electromagnetic fields / static charging (choose different location/switch off interfering device if possible)

Should other error messages occur, switch balance off and then on again. If the error message remains inform manufacturer.