<b>D</b>	<b>Installationsanleitung Akku</b>	Seite 2
<b>GB</b>	<b>Installation Information Rechargeable battery</b>	Page 6
<b>F</b>	<b>Notice d'Installation Accu</b>	Page 10
<b>E</b>	<b>Manual de Instalación Acumulador</b>	Página 14
<b>I</b>	<b>Istruzioni per l'installazione Accumulatore</b>	Pagina 18

---


---

# **KERN FEJ-A01**

Version 1.0 03/2006

---

---

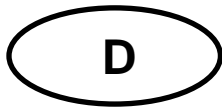# KERN FEJ-A01

Version 1.0 03/2006

## Installationsanleitung

### Akku

---

---

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---

#### 1 Technical data

- Operating duration approx. 6 h
- Charging time approx. 12 h via line adapter


#### 2 Scope of delivery

- Rechargeable battery
- 2 fastening screws
- Cover sheet with battery board


#### 3 To install rechargeable battery:

Remark: Simultaneous installation of the rechargeable battery pack and the relay exit is not possible.


- Disconnect the balance from the mains
- Remove weighing plate and carrier of weighing plate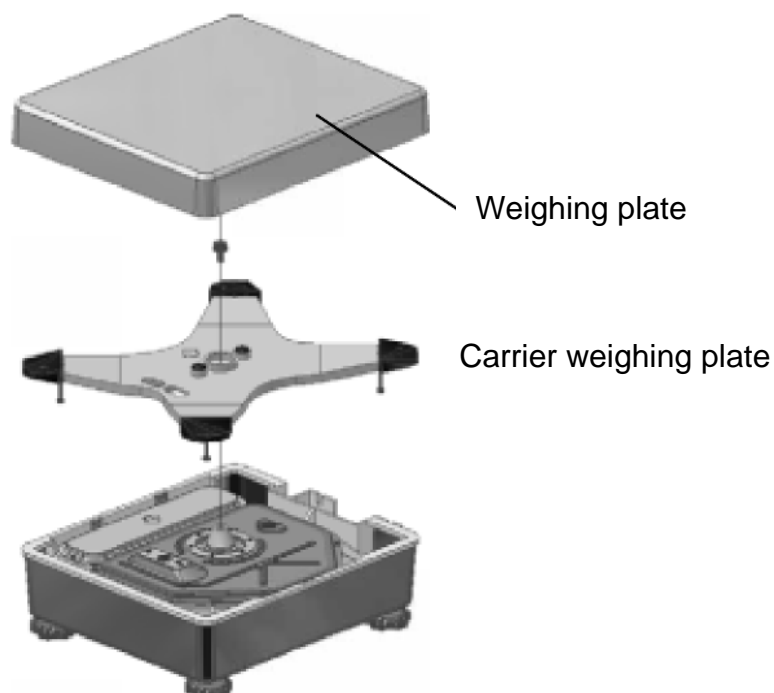

- Unscrew cover plate of battery compartment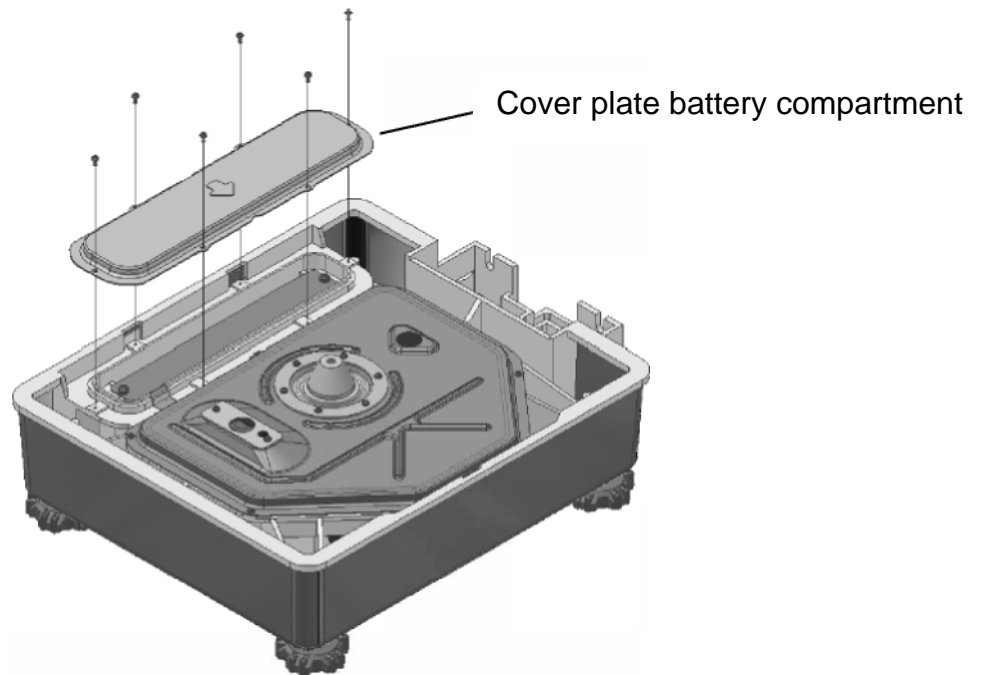


- Use screws to fix rechargeable battery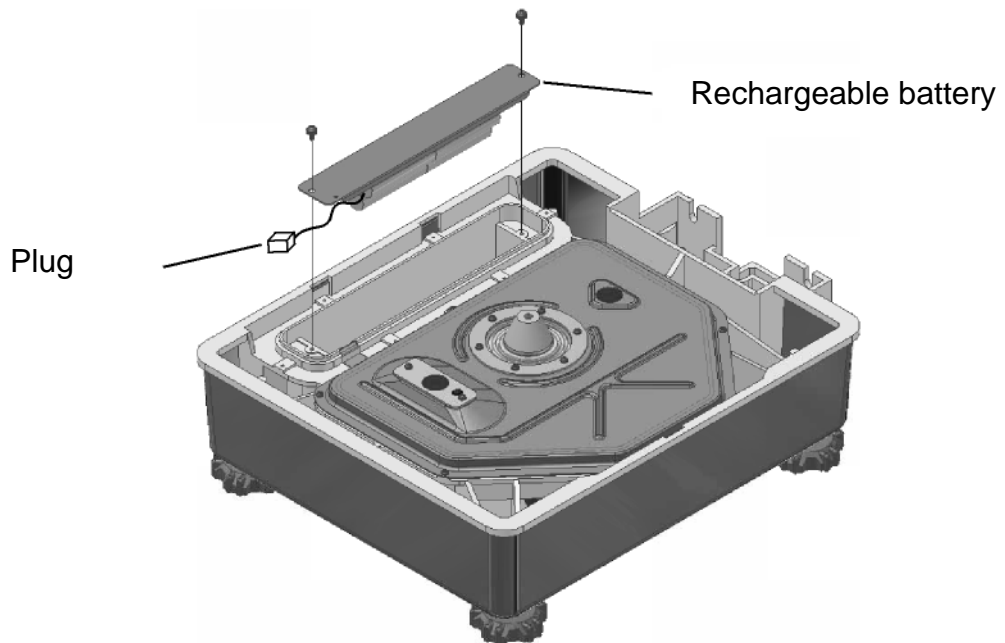

- Connect cable to plug
- Screw the cover plate of the battery compartment back on


#### 4 To install cover plate with battery board:

- Loosen nuts at rear end of display and remove front part of housing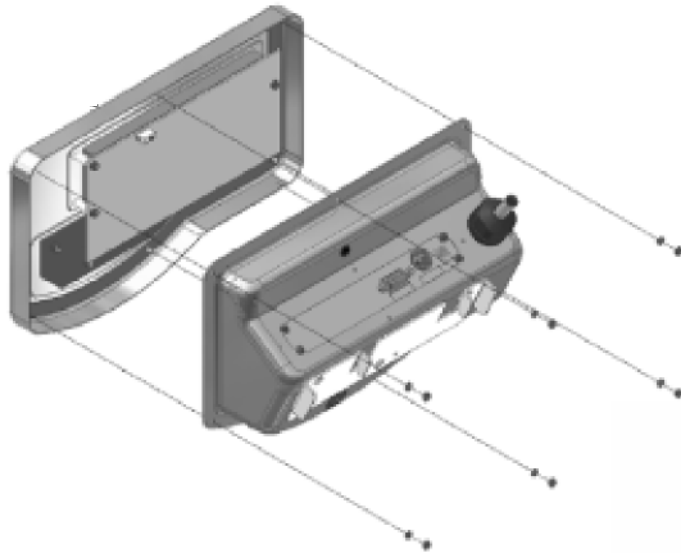

- Remove cover from rear part of housing
- Unscrew cover sheet with interface board and detach grounding cable connections CN 1 and CN2


Cover sheet with interface board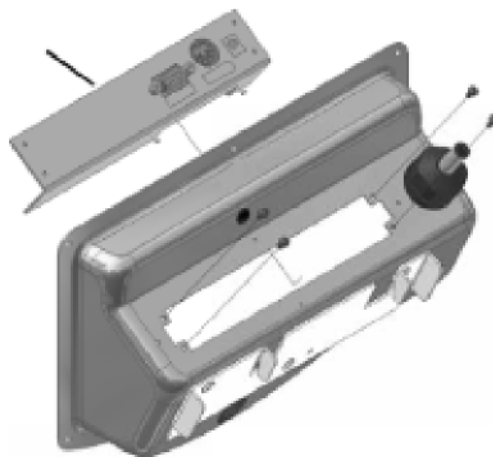

- Unscrew interface board from cover sheet and screw it to the cover sheet with battery board

- Screw down the cover sheet with battery board to the rear part of the housing

Cover sheet with  
battery board

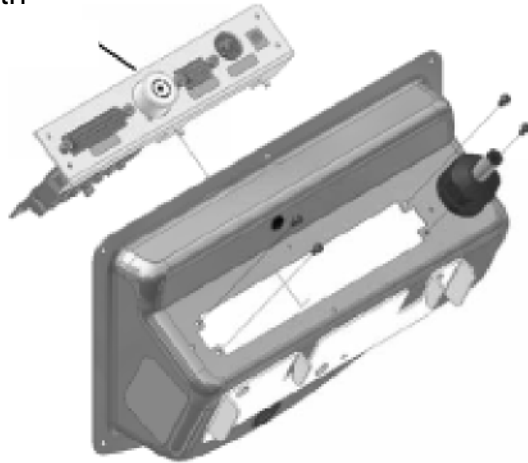- Plug in all connecting cables (CN 1, CN 2, CN 4) and screw the grounding cable back on
- Close housing

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