D

Ethernet-
schnittstelle

Seite 2

GB

Ethernet
interface

Page 6

F

Interface
Ethernet

Page 10

E

Interfaz
Ethernet

Página 14

I

Interfaccia
Ethernet

Pagina 18

KERN KMB-A01 / FTB-A09 / ITB-A17

Version 1.0 02/2007

ME Nr.: 22017111-

KMB-A01/ FTB-A09/ ITB-A17-IA-defsi-0710

KERN KMB-A01 / FTB-A09 / ITB-A17

Version 1.0 02/2007

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Version 1.0 02/2007

Ethernet interface

Supplement to operating instructions for balances with integrated Ethernet interface

If the balance has this built-in option, this additional interface may be used to couple the balance to the respective local network.

With the help of this interface the user can access the balance from the entire local network. As before, the same functionality (SICS commands) is available via the serial interface.

For details of these SICS interface commands please refer to the chapter "interface description" in the operating instructions for the balance.

For the connection to the network you will require the information and support of your local network administrator. The configuration shown below should be entered by your network administrator.

For this purpose, in the balance must be set the following adjustments.

The navigation through the menu can be seen in the balance instructions.

Installation:

Generally the Ethernet option under the menu item Option is shown as non-installable. This characterization applies to all options and is intended to provide the user with a quick overview of all available options.

COM 1		
COM 2		
COM 3		
Option	ETHERNET	not installed
	USB	not installed
	Analog	not installed

Notice: After the initial installation of the ETHERNET option, configuration is not possible before it has been enabled via the SICS command SOP.

The supply voltage for the Ethernet option, too, is turned on/off via the SICS command SOP (SET option). The syntax is shown in the overview below.

Command	Response	Description
SOP EO 1	SOP A	Enable Ethernet option
SOP EO 0	SOP A	Disable Ethernet option
SOP EO	SOP A 1 SOP A 0	Query setting Ethernet option enabled Ethernet option disabled

Serial interface:

The linking of the Ethernet option to the balance takes place internally via the serial interface COM 2 or the analogue option via COM 3. The picture below depicts the original settings. You may restore original settings by using COM Reset.

COM 1	Ethernet	
COM 2	Mode	Dialog
COM 3	Baud	9600
Option	Parity	8-none

Notice: The Ethernet option is supplied with the original setting 9600, 8-none. If the Ethernet option is installed or exchanged at a later date it is necessary to restore this original setting for the serial interface.

Net work:

Go to the menu item **Communication – Option – Ethernet** to set the parameters IP-address, subnet mask and gateway required for communication in the network. The following picture shows the original settings. Original settings may be restored by resetting the Ethernet.

Ethernet		
USB	IP-address	192.168.1.1
Analog	Subnet mask	255.255.255.0
	Gateway	0. 0. 0. 0

Attention: The quoted addresses are merely examples. You must ask your network administrator which settings are to be used.

Notice: “---.---.---.--- “ is displayed for the setting of the IP-address, subnet mask and gateway if no parameters are made available by the Ethernet option. Check the setting of the serial interface (original setting).

Check the correct installation of the Ethernet option in the balance.

IP-address:

The unique identification of a participant in the network is indicated by the IP-address. The IP-address is allocated to the participant either in a static or a dynamic way via DHCP.

DHCP is not supported by the Ethernet option. An IP-address consists of 4 bytes and is usually in decimal form. The byte value range is between 0 and 255. To achieve better readability, the bytes are separated by full stops.

Subnet Mask:

The IP-address comprises a Net-ID and a Host-ID. The Net-ID serves to address the network, the Host-ID to address the participant. Out of the 32 bits of the subnet mask the bits of the Net-ID are set to 1. Which part of the IP-address belongs to the Net-ID depends on the size of the network. The overview below depicts the 3 essential network categories.

Class	31	24	23	16	15	8	7	0
A	Net-ID			Host-ID				
B	Net-ID				Host-ID			
C	Net-ID						Host-ID	

It is possible to break up a network into further sub-networks (subnets). A part of the Host-ID is used to achieve this sub-division. However, this course of action reduces the number of potential participants in the subnet. In the original setting the subnet mask is set for a category 3 network.

Gateway :

The Net-ID allows the Ethernet option to detect whether the recipient of the data is registered in the same subnet. If this part of the recipient's IP-address matches the Ethernet option, the data will be directly forwarded to the recipient. If there is a difference, the recipient is not registered in the same subnet. The data will be forwarded for further transfer to the configurative gateway.

Course of action for testing communication:

After configuration a first test of the Ethernet option can be carried out. For simple tests and diagnosis in networks apply **Ping**. Ping allows you to test, whether a specific participant exists in the network and whether he/she is approachable. To activate the **Ping** command, go to the DOS input request (command prompt). The overview below shows the command syntax and possible responses.

Command	ping 192.168.1.1
Response	Response 192.168.1.1: bytes= 32, time= 2ms, TTL=64
No response	Timeout of response

Network connection:

If your PC is already connected via a switch / hub to the network, no further settings for the network card in the PC are required.

For a direct connection between Ethernet option and PC via a **crosspatch cable** comply with the settings for the Internet Protocol TCP/IP.

You find these settings under Windows via **Start-System Control-Network Connections**

To select features for the network card, click the right button of your mouse. Under **Internet Protocol TCP/IP Features** set options as follows:

Use the following IP-addresses:	
IP-address:	192.168. 1 . 2
Subnet mask :	255.255.255. 0
Standard gateway :	. . .

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