D

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Version 2.0 03/06

## Operating Instruction

Density determination set for precision weighing and analytical balances KERN ARS/ARJ/PRS/PRJ

---

---

### Contents:

<b>1</b>	<b>INTRODUCTION .....</b>	<b>12</b>
1.1	SCOPE OF DELIVERY .....	12
<b>2</b>	<b>EXPLANATORY NOTES TO DENSITY DETERMINATION .....</b>	<b>13</b>
2.1	METHODS OF DETERMINATION.....	13
2.1	DENSITY INDEX CALCULATION.....	13
2.2	SELECTING THE DENSITY DETERMINATION APPLICATION.....	13
2.3	CONFIGURING DENSITY DETERMINATION .....	14
2.4	STARTING AND INITIALISING THE DENSITY PROGRAM .....	14
2.5	DENSITY OF A LIQUID "LIQUID MODE" (WITH DENSITY SET KERN AR-A01) .....	15
2.6	DENSITY OF A SOLID „FIRMLY IN AIR MODE“.....	16
2.7	DENSITY OF A POROUS SOLID "FIRM POROUS MODE" .....	17
2.8	DENSITY OF A SOLID "FIRMLY ON FLOOR MODE" .....	19

# 1 Introduction

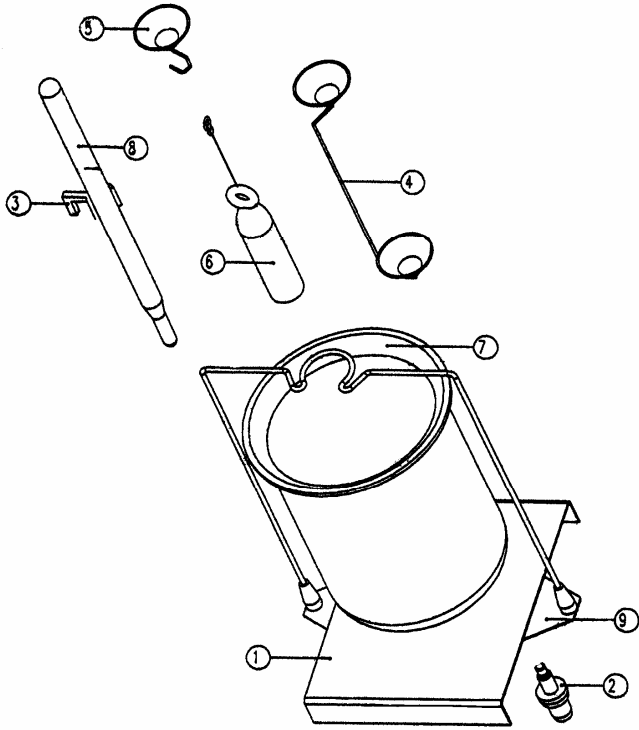
You can select between two different models when buying a density set as an accessory to your electronic balance:

**KERN AR-A01** For analytical balances of the **KERN ARS/ARJ** series

**KERN PR-A02** for analysing balances of the **KERN PRS/PLJ** series (only suitable for models with readability  $d = 0.001$  g).

These operating instructions only describe the operation of the density determination set. For further information on how to operate your balance please refer to the operating instructions supplied with each balance.

## 1.1 Scope of delivery

KERN AR-A01	KERN PR-A02												
 <ol style="list-style-type: none"> <li>1 Panel</li> <li>2 Cone for KERN ARS/ARJ density set (short cone)</li> <li>3 Support</li> <li>4 Twin beaker set</li> <li>5 Hook set</li> <li>6 Sinkers</li> <li>7 Glass beaker</li> <li>8 Thermometer</li> <li>9 Stand assembly</li> </ol>	 <table border="0"> <tr> <td>Platform</td> <td>KERN no. 904-101</td> </tr> <tr> <td>Beaker</td> <td>KERN no. 152-553</td> </tr> <tr> <td>Thermometer</td> <td>KERN no. 152-536</td> </tr> <tr> <td>Sinker</td> <td>KERN no. 152-532</td> </tr> <tr> <td>Filter bowl</td> <td>KERN no. 822-851</td> </tr> <tr> <td>Flush-mounted hook</td> <td>KERN no. 822-804</td> </tr> </table>	Platform	KERN no. 904-101	Beaker	KERN no. 152-553	Thermometer	KERN no. 152-536	Sinker	KERN no. 152-532	Filter bowl	KERN no. 822-851	Flush-mounted hook	KERN no. 822-804
Platform	KERN no. 904-101												
Beaker	KERN no. 152-553												
Thermometer	KERN no. 152-536												
Sinker	KERN no. 152-532												
Filter bowl	KERN no. 822-851												
Flush-mounted hook	KERN no. 822-804												

## 2 Explanatory notes to density determination

### 2.1 Methods of determination

To determine the density of solids and liquids, use the “density” application. There are various methods of determination to choose from:

- “LIQUID MODE“ Density of a liquid (density set AR-A01 only)
- “FIRMLY IN AIR MODE“ Density of a solid
- “FIRM POROUS MODE“ Density of a porous solid
- “FIRMLY ON FLOOR MODE“ Density of a solid with vessel on weighing tray

### 2.1 Density index calculation

The density index shows the percentage difference between two densities:

$$Index = \frac{Density\_1 \cdot Density\_2}{Density\_1} \cdot 100\%$$

To achieve this, the two densities are used in such a way that density<sub>1</sub> >= density<sub>2</sub>.

Either two successive densities are compared or the last density to be determined is compared to the reference density entered.

### 2.2 Selecting the density determination application

To activate the application menu press the «**MENU**» KEY and select „DENSITY“ application.

• SELECT APPLICATION		
SET APP:	OFF	Normal weighing mode
	...	...
	...	...
	DENSITY	Density determination
	...	...
	...	...

The sub-menus required to define density determination are now shown in the “SET APPLICATION“ menu.

## 2.3 Configuring density determination

• SETUP APPLICATION			
DENSITY	MODE	FIRMLY ON FLOOR	Solid and vessel on weighing tray
	MODE	FIRMLY IN AIR	Solid
	MODE	LIQUID	Liquids
	MODE	FIRM POROUS	Porous solids
	INDEX	ON/OFF	Index calculation ON/OFF
	REFERENCE	8.000	Reference density for INDEX calculation in g/cc (applied only if REFERENCE <> 0,000)
	TIME BASE	2,0	Time base for repetition in seconds (applied only if TIME BASE <> 0.0)
	REF. DENSITY	0,998205	Density of liquid used for measuring (default setting water at 20 °C)
TEMPERATURE	20 C	Temperature of water used for measuring in °C (REF. DENSITY is calculated accordingly)	

## 2.4 Starting and initialising the density program

To go to density program briefly press «».

If in "LIQUID MODE" this takes you directly to Density Determination. In all other modes the following message appears:

+	0,998205	g/cc	Current density of measured liquid
OK	KAL	T-H <sub>2</sub> O	20,0C Functions

### Occupancy of function keys:

Key	Functions
«OK»	Import density of measured liquid ==> Continue with relevant density determination.
«KAL»	Determine density of measured liquid. (See chapter 2.5. Density of liquid "liquid mode")
« T-H <sub>2</sub> O 20.0C »	Set reference density of water (currently set at: 20.0°C)
« T-H <sub>2</sub> O ---C »	Set reference density of water (currently not defined)

## 2.5 Density of a liquid “liquid mode“ (with density set KERN AR-A01)

This method determines the density of a liquid. Use a glass body of 10 cm<sup>3</sup> or 100 cm<sup>3</sup> volume.

### Display

### Key

### Step

	0,000	g
Air		SET

«T» Taring

Hook glass body into flush-mounted hook (fig. 1).

+	24,971	g
Air		SET

«⇔» Once value has stabilised, it is stored and displayed for 2 seconds

Place vessel with liquid on balance and submerge glass body without the glass body touching the bottom (fig. 2).

+	14,985	g
LIQUID		SET

«⇔» Once value has stabilised, it is stored and displayed for 2 seconds

+	0,9999	g/cc
DENSITY	<->	

Display of density of a liquid (as long as within 0.5 – 2.0 g/cc range).

+	0,20	%
INDEX	<->	

«↑» Display density index (only if INDEX ON).

	0,000	g
Air		SET

«T» Tare, balance switches to display in gram and is ready for the next density determination.

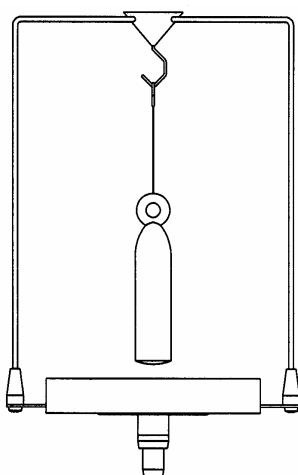
fig. 1

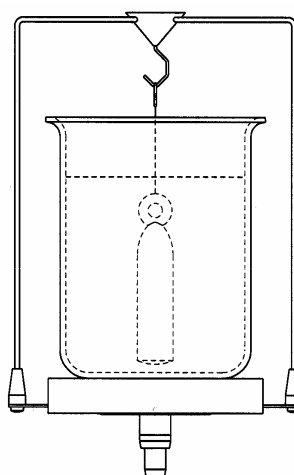
fig. 2

## 2.6 Density of a solid „firmly in air mode“

This method determines the density of a solid. The thermometer monitors the temperature of the measuring liquid in the vessel.

Display	Key	Step						
<table border="1"> <tr> <td>Air</td> <td>0,000</td> <td>g</td> </tr> <tr> <td></td> <td></td> <td>SET</td> </tr> </table>	Air	0,000	g			SET	«T»	Taring
Air	0,000	g						
		SET						
Put solid in top beaker (fig. 3 or fig. 3a).								
<table border="1"> <tr> <td>+</td> <td>24,972</td> <td>g</td> </tr> <tr> <td>Air</td> <td></td> <td>SET</td> </tr> </table>	+	24,972	g	Air		SET	«⇔»	Once value has stabilised, it is stored and displayed for 2 seconds
+	24,972	g						
Air		SET						
Put solid in bottom beaker (fig. 4 or fig. 4a).								
<table border="1"> <tr> <td>+</td> <td>16,982</td> <td>g</td> </tr> <tr> <td>LIQUID</td> <td></td> <td>SET</td> </tr> </table>	+	16,982	g	LIQUID		SET	«⇔»	Once value has stabilised, it is stored and displayed for 2 seconds
+	16,982	g						
LIQUID		SET						
<table border="1"> <tr> <td>+</td> <td>3,120</td> <td>g/cc</td> </tr> <tr> <td>DENSITY</td> <td>&lt;-&gt;</td> <td></td> </tr> </table>	+	3,120	g/cc	DENSITY	<->			Display of density of solid (as long as within 0.1 – 25.0 g/cc range).
+	3,120	g/cc						
DENSITY	<->							
<table border="1"> <tr> <td>+</td> <td>2,05</td> <td>%</td> </tr> <tr> <td>INDEX</td> <td>&lt;-&gt;</td> <td></td> </tr> </table>	+	2,05	%	INDEX	<->		«↑»	Display density index (only if INDEX ON).
+	2,05	%						
INDEX	<->							
<table border="1"> <tr> <td>Air</td> <td>0,000</td> <td>g</td> </tr> <tr> <td></td> <td></td> <td>SET</td> </tr> </table>	Air	0,000	g			SET	«T»	Tare, balance switches to display in gram and is ready for the next density determination.
Air	0,000	g						
		SET						

### Density determination by means of density set KERN AR-A01:

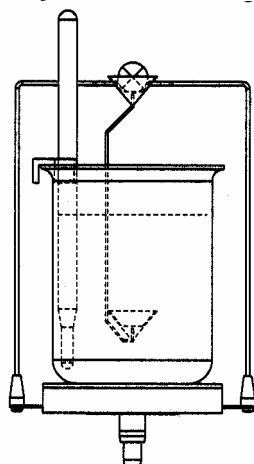
fig. 3

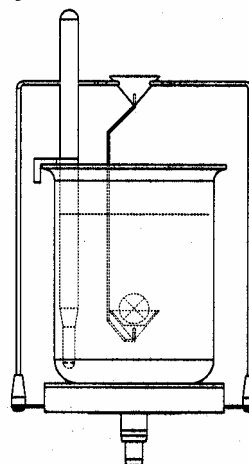
fig. 4

## Density determination by means of density set KERN PR-A02:

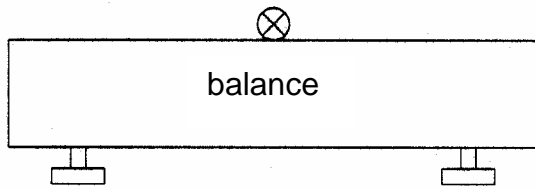
fig. 3a

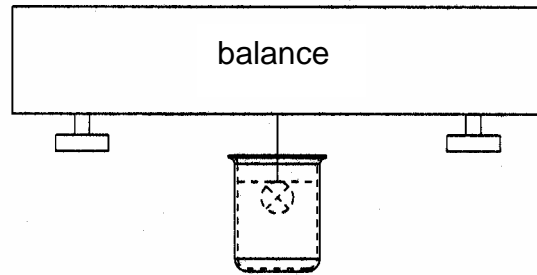
fig. 4a

### 2.7 Density of a porous solid "firm porous mode"

This method determines the density of a porous solid. The thermometer monitors the temperature of the measuring liquid in the vessel.

#### Display

#### Key

#### Step

	0,000	g
Air		SET

«T» Taring

Put porous solid in top beaker (fig. 5 or fig. 5a).

+	24,974	g
Air		SET

«⇔» Once value has stabilised, it is stored and displayed for 2 seconds

Seal pores of body, e. g. with wax. Now put solid in top beaker (fig. 6 or fig. 6a).

+	26,971	g
PORES CLOSED		SET

«⇔» Once value has stabilised, it is stored and displayed for 2 seconds

Put solid in bottom beaker (fig. 7 or fig. 7a)

+	16,984	g
LIQUID		SET

«⇔» Once value has stabilised, it is stored and displayed for 2 seconds

+	2,496	g/cc
DENSITY	<->	

Display of density of solid (as long as within 0.1 – 25.0 g/cc range).

+	2,05	%
INDEX	<->	

«↑» Display density index (only if INDEX ON).

	0,000	g
Air		SET

«T» Tare, balance switches to display in gram and is ready for the next density determination.

Density determination by means of density set KERN AR-A01:

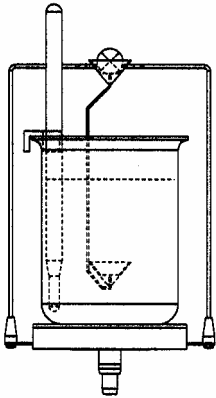
fig. 5

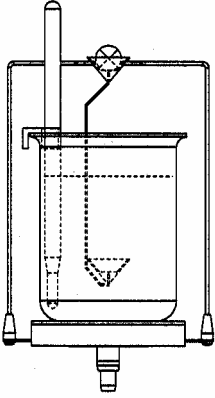
fig. 6

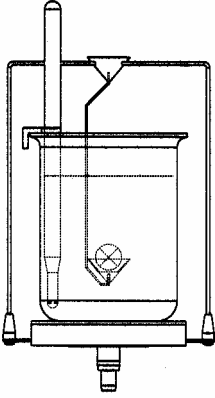
fig. 7

Density determination by means of density set KERN PR-A02:

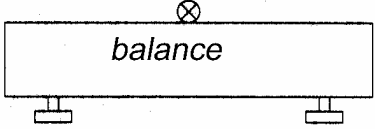
fig. 5a

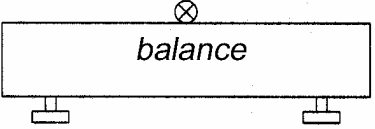
fig. 6a

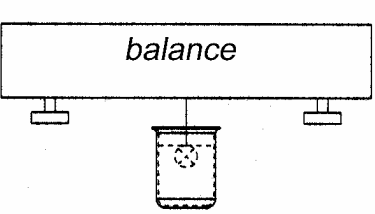
fig. 7a

## 2.8 Density of a solid “firmly on floor mode”

This method, too, determines the density of a solid; this time without the need for any contrivances. The thermometer monitors the temperature of the measuring liquid.

### Display

	0,000	g
FLOOR		SET

### Key

«T»

### Step

Put the temperate reference fluid (water) in a vessel, place it on the weighing platform and tare (fig. 8).

Put solid in vessel; positioned on floor (fig. 9).

+	24,971	g
FLOOR		SET

«⇨»

Once value has stabilised, it is stored and displayed for 2 seconds

Lift solid until it is no longer on the floor but remains completely submerged in the measuring fluid (Fig.10).

+	9,988	g
LIQUID		SET

«⇨»

Once value has stabilised, it is stored and displayed for 2 seconds

+	2,496	g/cc
DENSITY	<->	

Display of density of solid (as long as within 0.1 – 25.0 g/cc range).

+	7,36	%
INDEX	<->	

«↑»

Display density index (only if INDEX ON).

	0,000	g
FLOOR		SET

«T»

Tare, balance switches to display in gram and is ready for the next density determination.

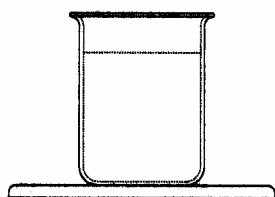
fig. 8

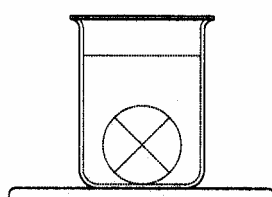
fig. 9

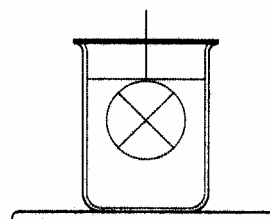
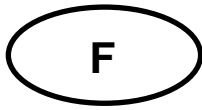
fig. 10

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