<b>D</b>	<b>Installationsanleitung Relais-Ausgang</b>	Seite 2
<b>GB</b>	<b>Installation Information Relay exit</b>	Page 9
<b>F</b>	<b>Notice d'Installation Sortie relais</b>	Page 16
<b>E</b>	<b>Manual de Instalación Salida de relais</b>	Página 23
<b>I</b>	<b>Istruzioni per l'installazione Uscitá su relé</b>	Pagina 30

---

---

# **KERN FEJ-A07**

Version 1.0 03/2006

---

---

# KERN FEJ-A07

Version 1.0 03/2006

## Installationsanleitung

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---

---

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---


#### Contents

<b>1</b>	<b><i>Installation</i></b>	<b>10</b>
<b>2</b>	<b><i>Specification</i></b>	<b>12</b>
2.1	Pin allocation of balance output plug D-SUB25P	12
2.2	Relay contact	13
2.3	Connection methods	13
2.4	Connection optoelectronic coupler	13
2.5	Relay input (example)	13
<b>3</b>	<b><i>Required menu settings</i></b>	<b>14</b>
<b>4</b>	<b><i>Operation</i></b>	<b>15</b>

## 1 Installation


Remark: Simultaneous installation of the rechargeable battery pack and the relay exit is not possible.

- Disconnect the balance from the mains!
- Loosen nuts at rear end of display and remove front part of housing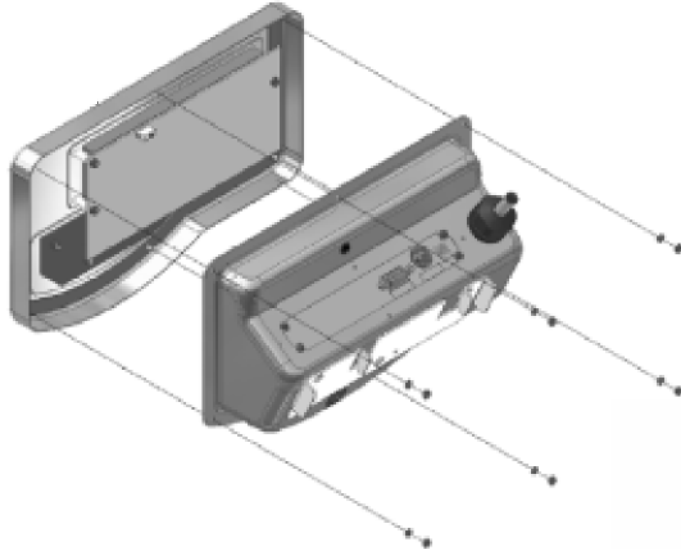

- Remove cover from rear part of housing
- Unscrew cover sheet with interface board and detach grounding cable connections CN 1 and CN2


Cover sheet with  
interface board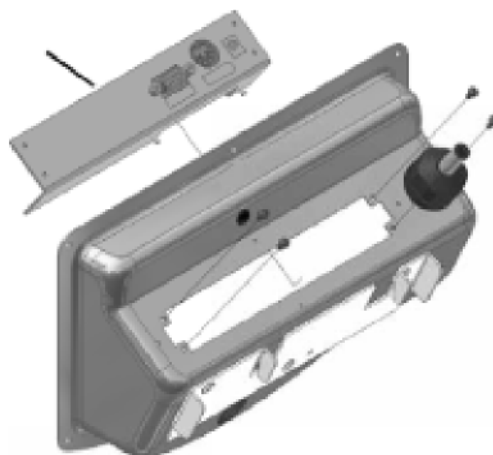

- Unscrew interface board from cover sheet and screw it to the relay cover sheet.

- Screw down the relay cover sheet to the rear part of the housing


Relay cover sheet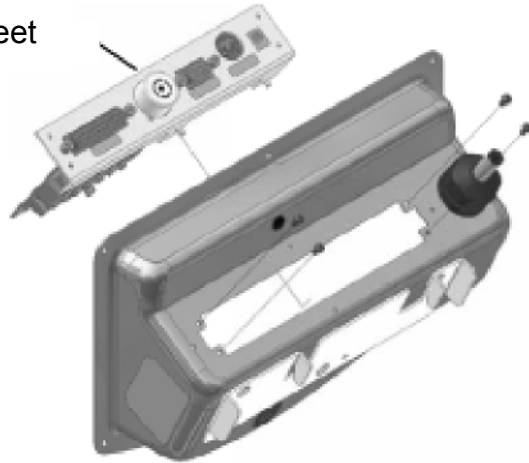

- Plug in all connecting cables (CN 1, CN 2, CN 4) and screw the grounding cable back on
- Close housing


## 2 Specification

### 2.1 Pin allocation of balance output plug D-SUB25P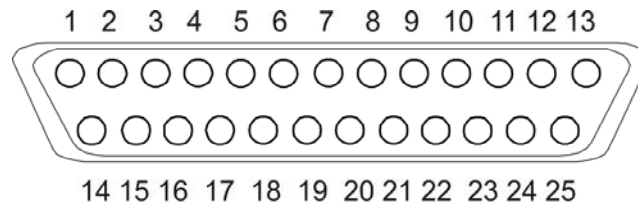

PIN	Signal		Input/ Output	Function
1	1. Limit (LOW)	NO	Output	ON, when first limit (LOW) is reached
2	1. Limit (LOW)	NC	Output	OFF, when first limit (LOW) is reached
3	1. Limit (LOW)	COM	-	Joint contact first limit (LOW)
4	2. Limit (OK)	NO	Output	ON, when second limit (OK) is reached
5	2. Limit (OK)	NC	Output	OFF, when second limit (OK) is reached
6	2. Limit (OK)	COM	-	Joint contact second limit (OK)
7	3. Limit (HIGH)	NO	Output	ON, when third limit (HIGH) is reached
8	3. Limit (HIGH)	NC	Output	OFF, when third limit (HIGH) is reached
9	3. Limit (HIGH)	COM	-	Joint contact third limit (HIGH)
10	4. Limit	NO	Output	ON, when forth limit is reached
11	4. Limit	NC	Output	OFF, when forth limit is reached
12	4. Limit	COM	-	Joint contact forth limit
13	5. Limit	NO	Output	ON, when fifth limit is reached
14	5. Limit	NC	Output	OFF, when fifth limit is reached
15	5. Limit	COM	-	Joint contact fifth limit
16	ERR	NO	Output	ON in case of error message
17	ERR	NC	Output	OFF in case of error message
18	ERR	COM	-	ERR joint contact
19	During measure- ment	NO	Output	ON, if measured quantity is 5 times greater than reading range
20	During measure- ment	NC	Output	OFF, if measured quantity is 5 times greater than reading range
21	During measure- ment	COM	-	Joint contact during measurement
22	Audio signal	NO	Output	On for audio signal
23	Audio signal	COM	-	Joint contact when audio signal is given
24	External signal (+)		Input	Contact for output control (+)
25	External signal (-)		Input	Contact for output control (-)

## 2.2 Relay contact

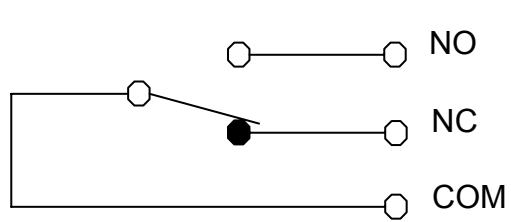
### Technical data:


AC 125 V/0.5A

DC 30V/1A


## 2.3 Connection methods

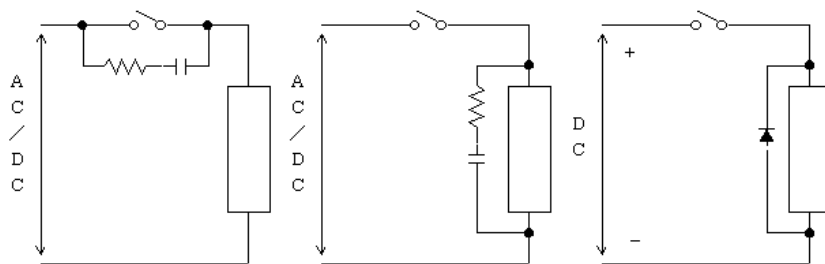## 2.4 Connection optoelectronic coupler
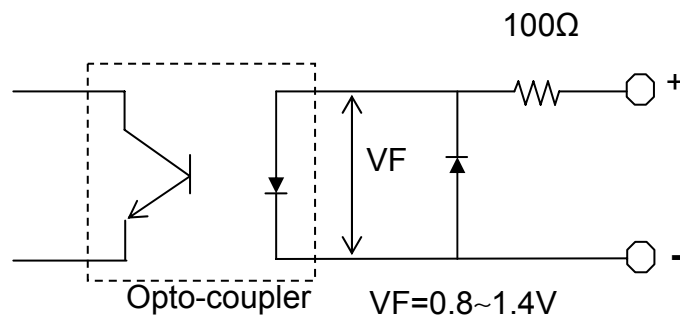
## 2.5 Relay input (example)
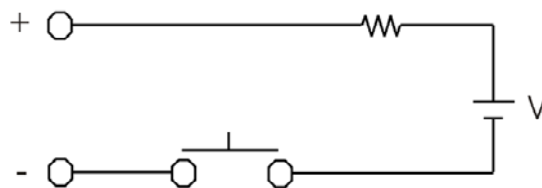
### 3 Required menu settings

For further information on menu settings please refer to the operating instructions supplied with all balances.

Default settings:

Additional functions	[2. SEL 2]	Tolerance weighing active
Display conditions of the tolerance marker	[21. Co. 1]	Tolerance marker is always displayed, even if standstill control is not yet displayed.
Tolerance range	[22. Li. 1]	Tolerance marker is displayed for the whole range.
Number of limiting points	[23. P1 2]	2- Limiting points (+/OK/-)
Assessment	[24. tYP. 1]	Evaluation for absolute values
Signal at limit 1	[25. bu1 0]	No signal at limit 1(-)
Signal at limit 2	[26. bu2 0]	No signal at limit 2(Ok)
Signal at limit 3	[27. bu3 0]	No signal at limit 3(+)
Signal at limit 4	[28. bu4 0]	No signal at limit 4
Signal at limit 5	[29. bu5 0]	No signal at limit 5
Display of Results	[2A. LG 1]	Display via +, OK or -
Output setting	[2b. r.o.c. 1]	Continuous output

## 4 Operation

- Connect mains adapter and start the balance by pressing the ON/OFF key
- Entering limiting values
- Evaluation of readings via relay contacts
- Status balance/contact

Status balance	Display	Contact
Reading outside condition	No reading display	inactive
Reading outside range	No reading display	inactive
Input of invalid values	Three ◀ symbols appear on the left	inactive
Error message (o-ERR/u-ERR)	o-ERR/u-ERR	<ul style="list-style-type: none"> <li>• OFF during measurement of limits 1-5</li> <li>• ON during ERR</li> </ul>
During input	Input appears on display	inactive
Waiting for stability message as per tares	[M ] flashing	Last display fixated
Stand-by	Stand by LED glowing	inactive
Sleep mode	Sleep LED glowing	Output of evaluation result
Automatic check of balance at starting	[00000000] glowing	inactive

- Output by external signal
  - ⇒ Contacts and audio signals are started by the weighing result, but may also be triggered by an external signal
  - ⇒ Whilst setting the menu [2b. r.o.c. 2] there is an output of the weighing result for approx. 200 ms after the signal was applied
  - ⇒ If a pulsed external signal is used, the output signal will follow after 100 ms at the earliest

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